00-8030 15 mL  
Histomount Mounting Solution  
Lot No.

00-8030 15 mL  
Histomount Mounting Solution  
Lot No.

### Histomount

### Histomount

#### INTRODUCTION

This product is an organic based mounting medium for immunohistochemical procedures. Organic insoluble chromogens such as DAB and BCIP/NBT are best preserved in this mounting medium.

#### REAGENT PREPARATION

This reagent is ready-to-use.

#### STORAGE

Store at room temperature away from flame.

#### RECOMMENDED GUIDELINES FOR USE

1. After immunostaining, dehydrate tissue sections in a graded series of alcohol and clear in xylene.
2. Cover tissue section with a few (2-4) drops of this mounting medium and put cover slip on.
3. Cap immediately after use.

Authorized Representative for IVDD 98/79/EC: Invitrogen Ltd.  
Inchinnan Business Park  
3 Fountain Drive  
Paisley  
PA4 9RF  
UK

#### INTRODUCTION

This product is an organic based mounting medium for immunohistochemical procedures. Organic insoluble chromogens such as DAB and BCIP/NBT are best preserved in this mounting medium.

#### REAGENT PREPARATION

This reagent is ready-to-use.

#### STORAGE

Store at room temperature away from flame.

#### RECOMMENDED GUIDELINES FOR USE

1. After immunostaining, dehydrate tissue sections in a graded series of alcohol and clear in xylene.
2. Cover tissue section with a few (2-4) drops of this mounting medium and put cover slip on.
3. Cap immediately after use.

Authorized Representative for IVDD 98/79/EC: Invitrogen Ltd.  
Inchinnan Business Park  
3 Fountain Drive  
Paisley  
PA4 9RF  
UK

For In Vitro Diagnostic Use. CAUTION: Not for human or animal therapeutic use.

For In Vitro Diagnostic Use. CAUTION: Not for human or animal therapeutic use.

[www.invitrogen.com](http://www.invitrogen.com)  
Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288  
E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)

[www.invitrogen.com](http://www.invitrogen.com)  
Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288  
E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)

PI008030

(Rev 06/09) DCC-09-0748

© 2007 Invitrogen Corporation. All right reserved. These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.

PI008030

(Rev 06/09) DCC-09-0748

© 2007 Invitrogen Corporation. All right reserved. These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.

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For In Vitro Diagnostic Use. CAUTION: Not for human or animal therapeutic use.

[www.invitrogen.com](http://www.invitrogen.com)  
Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288  
E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)

PI008030

(Rev 06/09) DCC-09-0748

Invitrogen Corporation. All right reserved. These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.

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Bevollmächtigter Repräsentant für IVDD 98/79/EC: Invitrogen Ltd.