
# Catalog Number: 195043, 1672346, 1672348 **Amphotericin B**

## Structure:

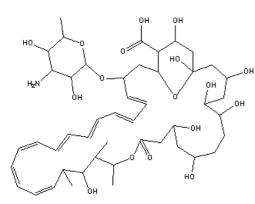

Molecular Formula: C<sub>47</sub>H<sub>73</sub>NO<sub>17</sub>

Molecular Weight: 924.11

CAS #: 1397-89-3

Synonyms: Fungizone; Amphozone; Fungilin; Ampho-Moronal

Source: Streptomyces species.

Physical Description: Yellow to orange powder. Depending on pH, amphotericin in solution is yellow to dark yellow.

Description: An antifungal agent. This product contains no animal-derived components.

Pka values: 5.5 and 10.0.3

**Solubility:** Soluble in acidic water (pH 2) or basic water (pH 11) (about 0.1 mg/ml). Water solubility can be increased with the addition of deoxycholic acid, sodium salt (increasing solubility with increasing concentration of deoxycholic acid). Soluble in Dimethylformamide (DMF) (2 to 4 mg/ml), DMF + 1 M HCI (3:1) (60 to 80 mg/ml), DMSO (30 to 40 mg/ml), propylene glycol; slightly soluble in methanol and methyl alcohol; insoluble in water, anhydrous alcohol, ether, benzene, and toluene. Solutions are stable for long periods between pH 4 and 10.<sup>1</sup>

Cat# 16723 is a concentrated solution of amphotericin at neutral to slightly basic pH, which is slightly cloudy due to the low solubility of amphotericin in aqueous buffer. The concentrated solution is stable at 2-8<sup>o</sup>C for up to 3 weeks. When diluted to the working concentration (see below for recommended concentration), the amphotericin will completely go into solution. Diluted amphotericin is stable for up to 7 days in culture at 37<sup>o</sup>C. Store diluted solutions aliquoted at -20°C or below for up to approximately 6 months.

Sterilization of solutions should be by filtration through teflon membrane filters; Amphotericin B should not be gamma irradiated or autoclaved.

#### Formulation (for 16723):

Component	mg/liter	Mol. Wt.	Mol. (mM)
Amphotericin B	0.25	924.1	0.00027
Sodium Deoxycholate	250	414.6	0.60

**Description:** An antimicrobial/antifungal agent effect against yeast and other fungi at a recommended concentration of 2.5 ug/L (prepared solution use 10 ml/L). It is a mixture of polyenes. It interrupts cell membrane permeability by binding sterols resulting in the loss of low molecular weight compounds from the cell. Amphotericin B may be toxic to some insect cell types. It is inactive against bacteria, rickettsia and viruses. It is typically stable in media at 37°C in the dark for approximately 3 days.<sup>21</sup>

#### Availability:

Catalog Number	Description	Size
195043	Amphotericin B, powder	100 mg
		250 mg

		500 mg 1 g 5 g
	Amphotericin B Solution, 250 ug/ml; formulated with deoxycholate	20 ml
1672348	and prepared in water for increased solubility, pH = 7-8	50 ml

### **References:**

- 1. Merck Index, 12th Ed., No. 627.
- 2. Pandey, R.C. and Rinehart, K.L., J. Antibiot., v. 29, 1035 (1976).
- 3. Mechlinski, et al., *Tetrahedron Letters*, 3873 (1970).
- 4. Holz, R.W., Antibiotics (N.Y.), v. 5, 313 (1979).
- 5. Medoff, G., et al., Ann. Rev. Pharmacol. Toxicol., v. 23, 303 (1983).
- 6. Clarkes's Isolation and Identification of Drugs, 2nd Ed., Moffatt, M.C., et al. (eds.); Pharmaceutical Press, p. 351 (1988).
- 7. Martindale, The Extra Pharmacopoeia, 30th Ed.; Pharmaceutical Press, pp. 315-318 (1993).
- 8. The Pharmacological Basis of Therapeutics, 8th Ed., Goodman and Gilman (eds.); Pergamon Press, p. 1165 (1990).
- 9. Data for Biochemical Research, 3rd Ed.; Oxford Press, pp. 296-297 (1989).
- 10. Analytical Profiles of Drug Substances, 6th Ed., Florey, K. (ed.); Academic Press, pp. 1-42 (1977).
- 11. Advances in Lipid Research, v. 14, 127 (1976).
- 12. Wildfeuer, A., et al., "In vitro activity of voriconazole against yeasts, moulds and dermatophytes in comparison with fluconazole, amphotericin B and griseofulvin." *Arzneimittelforschung*, v. 47:11, 1257-1263 (1997).
- 13. Kretschmar, M., et al., "Effects of amphotericin B incorporated into liposomes and in lipid suspensions in the treatment of murine candidiasis." *Arzneimittelforschung*, v. 46:7, 711-715 (1996).
- 14. Villani, P., et al., "Clinical and pharmacokinetic evaluation of a new lipid-based delivery system of amphotericin B in AIDS patients." *Arzneimittelforschung*, **v. 46:4**, 445-449 (1996).
- 15. Espinel-Ingroff, A., "Clinical utility of in vitro antifungal susceptibility testing." Rev. Esp. Quimioter., v. 13:2, 161-166 (2000).
- 16. Borrelli, P., et al., "Liposomal amphotericin B as first line and secondary prophylactic treatment for visceral leishmaniasis in a patient infected with HIV." *Ann. Ital. Med. Int.*, **v. 15:2**, 169-171 (2000).
- 17. Eldem, T. and Arican-Cellat, N., "High-performance liquid chromatographic determination of amphotericin B in a liposomal pharmaceutical product and validation of the assay." *J. Chromatogr. Sci.*, **v. 38:8**, 338-344 (2000).
- 18. Hahn, R.C. and Hamdan, J.S., "Effects of amphotericin B and three azole derivatives on the lipids of yeast cells of Paracoccidioides brasiliensis." *Antimicrob. Agents Chemother.*, **v. 44:7**, 1997-2000 (2000).
- 19. Younsi, M., et al., "Amphotericin B resistance and membrane fluidity in Kluyveromyces lactis strains." *Antimicrob. Agents Chemother.*, v. 44:7, 1911-1916 (2000).
- 20. Zager, R.A., "Polyene antibiotics: relative degrees of in vitro cytotoxicity and potential effects on tubule phospholipid and ceramide content." *Am. J. Kidney Dis.*, **v. 36:2**, 238-249 (2000).
- 21. Perlman, D., "Use of Antibiotics in cell culture media." *Methods in Enzymology: Cell Culture*, Jakoby, W.B. and Pastan, I.H. (eds.), Academic Press: New York, p. 112 (1979).