

E-Gel® Imager Light/Adaptor Base

Cat. nos. 4466602, 4466603, 4466604

Rev. Date: 6 August 2011

Part no. 25-1072 MAN0004954


Directions for installation of the E-Gel® Safe Imager™, E-Gel® iBase™ Device, and E-Gel® Go! Base with the E-Gel® Imager UV/Blue-Light Base, or E-Gel® Imager Adaptor Base are described below. For detailed instructions refer to the **E-Gel® Imager Gel Documentation System** manual included in the software CD, or contact Technical Support.

E-Gel® Imager Light Base



ON/OFF
Power Button


E-Gel® Imager Light Base Ports

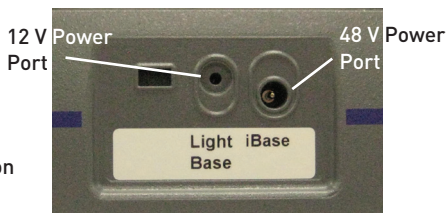

E-Gel® Imager Adaptor Base



E-Gel® Imager Adaptor Base Port

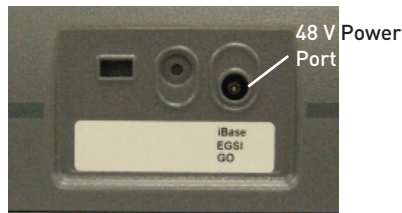

Kit contents

Catalog no.	Components	Quantity
4466602	E-Gel® Imager UV Light Base Universal 12 V Power Cord	1 1
4466603	E-Gel® Imager Blue-Light Base Universal 12 V Power Cord	1 1
4466604	E-Gel® Imager Adaptor Base DC Connector Cable	1 1

Intended Use

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

Installing the E-Gel® iBase™ with the E-Gel® Imager Light Base


The E-Gel® Imager Blue-Light Base and E-Gel® Imager UV Light Base can be used to provide transillumination for the E-Gel® iBase™.

For best results, affix the E-Gel® Imager Light Diffuser to the bottom of the E-Gel® iBase™ when using this configuration.



1. Place the E-Gel® iBase™ directly onto the transilluminator surface of the E-Gel® Imager Light Base.
2. Connect the DC Connector Cable between the E-Gel® iBase™ and the E-Gel® Imager Adaptor Base.

E-Gel® Imager Adaptor Base

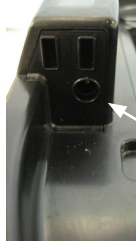

DC Connector Cable



E-Gel® iBase™ Device




3. Connect the E-Gel® iBase™ power supply cable to the 48 V power port at the rear of the E-Gel® Imager Light Base.
4. The E-Gel® iBase™ Device is now ready for use with the E-Gel® Imager system.

Installing the E-Gel® iBase™ with the E-Gel® Imager Adaptor Base

The E-Gel® Safe Imager™ can be used with the E-Gel® Imager Adaptor Base to provide transillumination for imaging with the E-Gel® iBase™ Device.

Note: Make sure to press the ON button of the E-Gel® Safe Imager™ for 2 seconds to extend the duration of transillumination.

Important: There are two types of plug on the E-Gel® Safe Imager™. Choose the appropriate installation method for the E-Gel® iBase™ / E-Gel® Safe Imager™ based on whether your E-Gel® Safe Imager™ has Plug A, or Plug B.

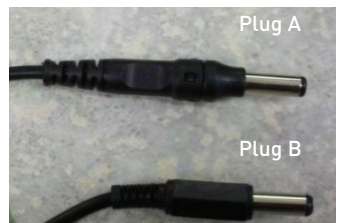

Plug A Installation Method

1. Place the E-Gel® Safe Imager™ into the well of the E-Gel® Imager Adaptor Base.
2. Connect the E-Gel® Safe Imager™ (**Plug A**) to the E-Gel® Imager Adaptor Base.
3. Connect the DC Connector Cable between the E-Gel® Safe Imager™ and the E-Gel® iBase™ Device.



E-Gel® Imager Adaptor Base



E-Gel® Safe Imager™



DC Connector Cable



E-Gel® iBase™ Device



4. Connect the E-Gel® iBase™ power supply cable to the 48 V power port at the rear of the E-Gel® Imager Adaptor Base.
5. The E-Gel® iBase™ Device is now ready for use with the E-Gel® Imager system.

Plug B Installation Method

1. Place the E-Gel® Safe Imager™ into the well of the E-Gel® Imager Adaptor Base.
2. Connect the DC Connector Cable between the E-Gel® Safe Imager™ and the E-Gel® Imager Adaptor Base.
3. Connect the E-Gel® Safe Imager™ (**Plug B**) to the E-Gel® iBase™ Device.

E-Gel® Imager Adaptor Base



DC Connector Cable



E-Gel® Safe Imager™



E-Gel® iBase™ Device



4. Connect the E-Gel® iBase™ power supply cable to the 48 V power port at the rear of the E-Gel® Imager Adaptor Base.
5. The E-Gel® iBase™ Device is now ready for use with the E-Gel® Imager system.

Installing the E-Gel® Go! Base with the E-Gel® Imager Adaptor Base


The E-Gel® Go! Base can be used with the E-Gel® Imager Adaptor Base either with or without the E-Gel® Go! Portable Battery Pack.

When not using the Portable Battery Pack, the E-Gel® Imager Adaptor Base can be used to provide power for the E-Gel® Go! Base.



1. Place the E-Gel® Go! Base into the well of the E-Gel® Imager Adaptor Base on the dedicated peg niches.
2. Connect the DC Connector Cable between the E-Gel® Go! Base and the E-Gel® Imager Adaptor Base.

E-Gel® Imager Adaptor Base

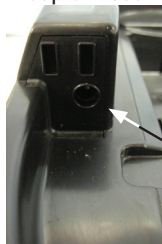

DC Connector Cable



E-Gel® iBase™ Device



3. Connect the E-Gel® Go! Base power supply cable to the 48 V power port at the rear of the E-Gel® Imager Adaptor Base.
4. The E-Gel® Go! Base is now ready for use with the E-Gel® Imager system.

Limited Use Label License: Research Use Only: The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial services of any kind, including, without limitation, reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies, 5791 Van Allen Way, Carlsbad, California 92008.

Limited Use Label License

©2011 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

For technical support, email tech_support@invitrogen.com.
For country-specific contact information, visit www.invitrogen.com.
To reorder visit www.invitrogen.com

