

## PRODUCT INFORMATION

# Thermo Scientific TurboFect *in vivo* Transfection Reagent

120 µl (for 20 transfections using 50 µg DNA)

# \_\_\_\_\_

Lot: \_\_\_\_\_

Expiry Date: \_\_\_\_\_

Store at 4°C



[www.thermoscientific.com/onebio](http://www.thermoscientific.com/onebio)

## Description

Thermo Scientific TurboFect *in vivo* Transfection Reagent is a sterile solution of a cationic polymer in water. The polymer forms compact, stable, positively charged complexes with DNA. These complexes protect DNA from degradation and facilitate gene delivery *in vivo*. Reagent causes no detectable inflammatory response and is suitable for DNA administration via various routes, including intravenous, intraperitoneal, intratumoral injection, etc.

## Reagents to be Supplied by the User

Sterile 5% w/v glucose solution for dilution of DNA.

## General Considerations

### DNA Quality Requirements

DNA quality is critical for successful transfection. An  $A_{260}/A_{280}$  ratio of 1.8 or higher is recommended. DNA should be sterile and free of any contaminants such as endotoxins.

### Choice of promoter

High gene expression depends on both the promoter under which the gene of interest is expressed and the targeted tissue/organ. Cytomegalovirus (CMV) promoter is best known for high gene expression in a wide variety of cells. Some researchers prefer LTR, simian virus (SV40) or Rous sarcoma virus (RSV) promoters.

### Transfection Reagent/DNA Ratio

The volume of transfection reagent used depends on the amount of DNA, transgene and animals to be transfected. The ratios presented in the protocols below are starting amounts and can be further optimized for best results.

### Amount of DNA and suggested injection volume

The amount of DNA and suggested injection volume depend on the experimental animal and the route of administration (see Table 1) as well as the targeted tissue or organ and the expression vector.

To prevent precipitation of TurboFect™/DNA complexes, the final concentration of DNA in the injection mix should not exceed 0.5 µg/µl.

## General Protocol for *in vivo* transfection

The protocol is optimized for transfection of mice via the tail vein.

Quantities and volumes should be scaled-up according to the animal to be transfected (see Table 1).

1. Dilute 50 µg of DNA in 400 µl sterile 5% glucose solution. Vortex gently and spin down briefly.
2. Briefly vortex TurboFect *in vivo* Transfection Reagent and add 6 µl of it to the diluted DNA. Mix immediately by pipetting or vortexing.
3. Incubate for 15-20 min. at room temperature.
4. Perform injections.
5. Monitor gene expression with the method most suitable for your studies.

### Note

The  $A_{260}/A_{280}$  ratio should be at least 1.8 for purified DNA. It is important to use endotoxin-free DNA (less than 0.1 EU/1 µg DNA).

The amount of DNA and suggested injection volume depend on the experimental animal and the route of administration (see Tables 1 and 2) as well as on the targeted tissue or organ and the expression vector.

(continued on reverse page)

**Table 1. Suggested amount of DNA and injection volume**

Animal	Route of injection	Suggested amount of DNA ( $\mu\text{g}$ )	Suggested injection volume
Adult Mouse	Intravenous (tail)	50	400 $\mu\text{l}$
	Intraperitoneal	100	800 $\mu\text{l}$
	Intratumoral	10	50 $\mu\text{l}$
	Brain	2.5	5 $\mu\text{l}$
Newborn mouse	Brain	1	2 $\mu\text{l}$
Nude mouse	Intravenous (tail)	50	400 $\mu\text{l}$
	Intratumoral	10	50 $\mu\text{l}$
Adult rat	Intravenous (tail)	500	2 ml
	Brain	20	25 $\mu\text{l}$
Adult rabbit	Intravenous	5000	10 ml

**Table 2. Scale-up ratios for transfection using TurboFect *in vivo* Transfection Reagent**

Amount of DNA ( $\mu\text{g}$ )	Volume of TurboFect <i>in vivo</i> Transfection Reagent ( $\mu\text{l}$ )	
	Recommended	Range
1	0.12	0.1-0.16
5	0.6	0.5-0.8
10	1.2	1-1.6
50	6	5-8

**CERTIFICATE OF ANALYSIS**

Transfection efficiency was tested on BALB/c female mice (7-8 weeks old) using 50  $\mu\text{g}$  Firefly luciferase coding plasmid and 6  $\mu\text{l}$  of TurboFect *in vivo* Transfection reagents per mouse. The delivery of the transgene was evaluated by measurement of Firefly luciferase activity in total protein extracted from the lungs of transfected animals. 100% of tested mice expressed the Firefly luciferase.

Quality authorized by:

 Jurgita Zilinskiene

## Troubleshooting

Problem	Possible Cause and Solution
<b>Low transfection efficiency</b>	<b>Suboptimal reagent/DNA ratio.</b> Optimize the amount of transfection reagent added to the fixed amount of DNA. <b>Suboptimal quantity of DNA.</b> Optimize the amount of DNA used for transfection. Keep the transfection reagent/DNA ratio constant. <b>Poor DNA quality.</b> Use high quality endotoxin-free DNA with an $A_{260}/A_{280}$ ratio greater than 1.8.
<b>Mortality</b>	<b>Suboptimal reagent/DNA ratio.</b> Decrease the amount of transfection reagent added to the fixed amount of DNA. <b>Suboptimal quantity of DNA.</b> Decrease the amount of DNA. Keep the transfection reagent/DNA ratio constant. <b>Poor DNA quality.</b> Use high quality endotoxin-free DNA with an $A_{260}/A_{280}$ ratio greater than 1.8.

### Note

This product or the use of this product is covered by US patent application US20100041739A1 and corresponding counterparts. The purchase of this product includes a non-transferable license to use this product for the purchaser's internal research. All other commercial uses of this product, including without limitation product use for diagnostic purposes, resale of product in the original or any modified form or product use in providing commercial services require a separate license. For further information on obtaining licenses please contact [info.fermentas.lt@thermofisher.com](mailto:info.fermentas.lt@thermofisher.com)

### **PRODUCT USE LIMITATION**

This product is developed, designed and sold exclusively for research purposes and in vitro use only. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans.

Please refer to [www.thermoscientific.com/onebio](http://www.thermoscientific.com/onebio) for Material Safety Data Sheet of the product.

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.