

ELISA for Mouse IL-4

Product Code: 3311-1H-6

CONTENTS, development kit for 6 plates:

Vial 1 (red top)

Monoclonal antibody 11B11 (150 μ l)

Concentration: 1 mg/ml

Vial 2 (blue top)

Biotinylated monoclonal antibody BVD6-24G2 (80 μ l).

Concentration: 1 mg/ml

Vial 3 (white top)

Streptavidin-Horseradish Peroxidase (80 μ l)

Vial 4

Recombinant mouse IL-4 standard (1 μ g)

To ensure total recovery of stated quantity, vials have been overfilled.

STORAGE:

Shipped at ambient temperature. On arrival box 1 should be stored refrigerated at 4-8°C and box 2 should be stored at -20°C.

General

Intended use: For quantitative determination of native mouse IL-4 in solution, e.g. cell culture supernatant and serum/plasma samples.


Reagents: Antibodies are supplied in sterile-filtered (0.2 μm) PBS with sodium azide (0.02%). Streptavidin-HRP is supplied in PBS with 1% BSA and 0.15% Kathon CG.

Standard range: 4-400 pg/ml

Intra-assay variation: < 5%

Standard calibration: 1 ng of supplied standard equals 37 U of 91/656 NIBSC*-standard according to repeated calibrations. Calibration is batch-specific.

*National Institute of Biological Standards and Control, UK.

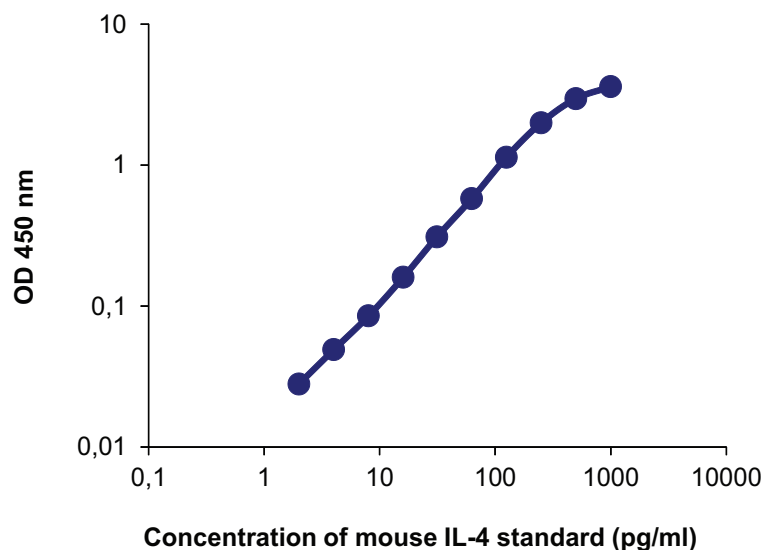

Guidelines for Mouse IL-4 ELISA

- Day 1**
1. Dilute mAb 11B11, diluted to 2 $\mu\text{g}/\text{ml}$ in PBS, pH 7.4, and filter the solution through a 0.2 μm filter. Coat a high protein binding ELISA plate with the solution by adding 100 $\mu\text{l}/\text{well}$. Incubate overnight at 4-8°C.
- Day 2**
2. Wash twice with PBS (200 $\mu\text{l}/\text{well}$).
 3. Block plate by adding 200 $\mu\text{l}/\text{well}$ of PBS with 0.05% Tween 20 containing 0.1% BSA (incubation buffer). Incubate for 1 hour at room temperature.
 4. Wash five times with PBS containing 0.05% Tween.
 5. Prepare mouse IL-4 standard by reconstituting contents of vial 4 in 1 ml PBS to make a stock solution of 1 $\mu\text{g}/\text{ml}$. Allow the standard to dissolve for 5 minutes, mix thoroughly and aliquot. Store at -20°C and avoid repeated freeze-thaw cycles of the standard aliquots. For the test, prepare dilutions of the stock using the standard range as a guideline.
 6. Add 100 $\mu\text{l}/\text{well}$ of samples or standards diluted in incubation buffer and incubate for 2 hours at room temperature.
 7. Wash as in step 4.
 8. Add 100 $\mu\text{l}/\text{well}$ of mAb BVD6-24G2-biotin at 0.1 $\mu\text{g}/\text{ml}$ in incubation buffer. Incubate for 1 hour at room temperature.
 9. Wash as in step 4.
 10. Add 100 $\mu\text{l}/\text{well}$ of Streptavidin-HRP diluted 1:1000 in incubation buffer. Incubate for 1 hour at room temperature. **Please note that sodium azide used in buffers will inhibit HRP activity.**
 11. Wash as in step 4.
 12. Add 100 $\mu\text{l}/\text{well}$ of appropriate substrate solution.
 13. Measure the optical density in an ELISA reader after suitable developing time. If required stop the reaction first.

NOTE; for research use only.

MABTECH shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.



MABTECH AB
Box 1233
SE-131 28 Nacka Strand
Sweden
Tel: +46 8 716 27 00
Fax: +46 8 716 27 01
E-mail: mabtech@mabtech.com
www.mabtech.com

MABTECH Inc
M.E.B. 220
3814 West Street
Cincinnati, OH 45227
USA
Tel: +1 513 871 4500
Fax: +1 513 871 7353
E-mail: mabtech.usa@mabtech.com

MABTECH AB Büro Deutschland
Germany
Tel: +49 40 4135 7935
Fax: +49 40 4135 7945
E-mail: mabtech.de@mabtech.com

2012-02-03

Developed and manufactured by MABTECH AB, Sweden, whose quality management system complies with the following standards:



MABTECH AUSTRALIA Pty Ltd
44 Gresswell Road
Macleod, VIC 3085
Australia
Tel: +61 3 9459 9630
Fax: +61 3 9455 0084
E-mail: mabtech.au@mabtech.com

MABTECH AB Bureau de liaison France
BP 255, 1300 route des Crêtes
06905 Sophia Antipolis
France
Tel: +33 (0)4 92 38 80 70
Fax: +33 (0)4 92 38 80 71
E-mail: mabtech.fr@mabtech.com