

TMB

Product code:	3651-10
Contents:	Contains tetramethylbenzidine, TMB, 1.25 mM in acetate buffer, pH 4.9. Also contains stabilizers and precipitating agents.
Applications:	Substrate for horseradish peroxidase conjugates in ELISpot assay.
Instructions:	Filter the ready-to-use substrate solution (TMB) through a 0.45 µm filter and add to ELISpot plates after incubation with horseradish peroxidase and subsequent washing. Develop until spots emerge. Stop the colour development by washing extensively in tap water. Use water of good quality since poor quality may cause fading of TMB spots.
Storage:	Shipped at ambient temperature. On arrival store substrate refrigerated at 4-8°C. Warm substrate to reaction temperature before use.



Note; for research use only.

Mabtech shall not be liable for the use or handling of the product or for consequential, special, indirect or incidental damages therefrom.



MABTECH AB
Box 1233
SE-131 28 Nacka Strand
Sweden
Tel +46 8 716 27 00
Fax +46 8 716 27 01
mabtech@mabtech.com
www.mabtech.com

Mabtech, Inc.
M.E.B. Suite 220
3814 West Street
Mariemont, OH 45227
USA
Toll Free: +1 866 ELI-SPOT
Tel: +1 513 871 4500
Fax: +1 513 871 7353
mabtech.usa@mabtech.com

MABTECH AUSTRALIA
Pty Ltd
resolvingIMAGES
Unit 22
Thomastown Victoria 3074
Australia
Tel: +61 3 9466 4007
Fax: +61 3 9466 4003
mabtech.au@mabtech.com

Mabtech AB
Büro Deutschland
GERMANY
Tel: +49 40 4135 7935
Fax: +49 40 4135 7945
mabtech.de@mabtech.com

Mabtech AB
Bureau de liaison France
BP 255, 1300 route des
Crêtes
06905 Sophia Antipolis
FRANCE
Tel: +33 4 92 38 80 70
Fax: +33 4 92 38 80 71
mabtech.fr@mabtech.com