

Technical Data Sheet

V500 Mouse Anti-Human CD3

Product Information


Material Number:	560772
Size:	25 tests
Vol. per Test:	5 µl
Clone:	SP34-2
Isotype:	Mouse IgG1, λ
Reactivity:	Human
	QC Testing: baboon or rhesus or cynomolgus
Storage Buffer:	Aqueous buffered solution containing protein stabilizer and ≤0.09% sodium azide.

Description

Clone SP34-2 is a mouse IgG1 isotype monoclonal antibody, descendant of SP34 (mouse IgG3), with the same specificity and reactivity pattern as the parent clone. It cross-reacts with a major subset of peripheral blood lymphocytes, but not monocytes or granulocytes, of baboon, and rhesus, cynomolgus, and pigtail macaque monkeys. The distribution on lymphocytes is similar to that observed with normal human donor lymphocytes with the majority of CD3-positive cells being negative when dual stained with antibodies to B or NK cells markers. SP34-2 is also capable of inducing cell proliferation on both human and non-human primate PBMC.

The antibody is conjugated to BD Horizon™ V500, which has been developed for use in multicolor flow cytometry experiments and is available exclusively from BD Biosciences. It is excited by the Violet laser with an Ex max of 415 nm and Em Max at 500 nm. BD Horizon V500 conjugates emit at a similar wavelength to Amcyan yet exhibit reduced spillover into the FITC channel. For more information on BD Horizon V500, visit bdbiosciences.com/colors.

When compensating dyes in this spectral range (such as Horizon™ V500 and AmCyan), the most accurate compensation can be obtained using single stained cellular controls. Due to spectral differences between cells and beads in this channel, using BD CompBeads can result in spillover errors for V500 and AmCyan reagents. Therefore, the use of BD CompBeads or BD CompBeads Plus to determine spillover values for these reagents is not recommended. Different V500 reagents (e.g. CD4 vs. CD45) can have slightly different fluorescence spillover therefore, it may also be necessary to use clone specific compensation controls when using these reagents.

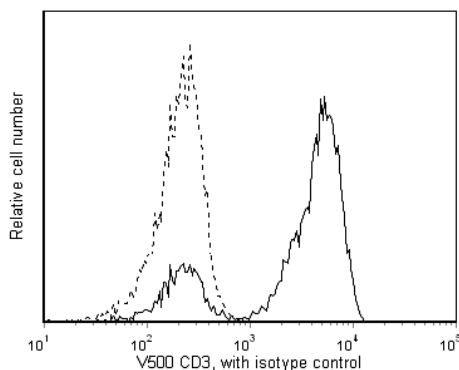

Analysis of CD3 on peripheral blood lymphocytes of rhesus macaque (*Macaca mulatta*) monkey. Peripheral blood lymphocytes were stained with BD Horizon™ V500 Mouse Anti-Human CD3 (solid line) and compared to whole blood stained with BD Horizon™ V500 Mouse IgG1, κ Isotype Control (clone X40, Cat. No. 560787, dotted line). Lymphocytes were selected by light scatter profile. Flow cytometry was performed on a BD FACSCanto™ II flow cytometry system.

Preparation and Storage

The monoclonal antibody was purified from tissue culture supernatant or ascites by affinity chromatography.

The antibody was conjugated with BD Horizon™ V500 under optimum conditions, and unreacted BD Horizon™ V500 was removed.

Store undiluted at 4°C and protected from prolonged exposure to light. Do not freeze.

BD Biosciences

bdbiosciences.com

United States	Canada	Europe	Japan	Asia Pacific	Latin America/Caribbean
877.232.8995	888.268.5430	32.53.720.550	0120.8555.90	65.6861.0633	0800.771.7157

For country-specific contact information, visit bdbiosciences.com/how_to_order/

Conditions: The information disclosed herein is not to be construed as a recommendation to use the above product in violation of any patents. BD Biosciences will not be held responsible for patent infringement or other violations that may occur with the use of our products. Purchase does not include or carry any right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of Becton Dickinson and Company is strictly prohibited.

For Research Use Only. Not for use in diagnostic or therapeutic procedures. Not for resale.

BD, BD Logo and all other trademarks are the property of Becton, Dickinson and Company. ©2011 BD



Application Notes

Application

Flow cytometry

Routinely Tested

Suggested Companion Products

Catalog Number	Name	Size	Clone
560787	V500 Mouse IgG1, κ Isotype Control	0.1 mg	X40

Product Notices

1. This reagent has been pre-diluted for use at the recommended Volume per Test. We typically use 1×10^6 cells in a 100- μ l experimental sample (a test).
2. Please refer to www.bdbiosciences.com/pharming/en/protocols for technical protocols.
3. BD Horizon™ V500 has a maximum absorption of 415 nm and maximum emission of 500 nm. Before staining with this reagent, please confirm that your flow cytometer is capable of exciting the fluorochrome and discriminating the resulting fluorescence.
4. Caution: Sodium azide yields highly toxic hydrazoic acid under acidic conditions. Dilute azide compounds in running water before discarding to avoid accumulation of potentially explosive deposits in plumbing.
5. For fluorochrome spectra and suitable instrument settings, please refer to our Fluorochrome Web Page at www.bdbiosciences.com/colors.

References

Bleavins MR, Brott DA, Alvey JD, de la Iglesia FA. Flow cytometric characterization of lymphocyte subpopulations in the cynomolgus monkey (*Macaca fascicularis*). *Vet Immunol Immunopathol.* 1993; 37(1):1-13. (Biology)

Conrad ML, Davis WC, Koop BF. TCR and CD3 antibody cross-reactivity in 44 species. *Cytometry A.* 2007; 71(11):925-933. (Biology)

Engel P, Wagner N, Zhou L, et al. CD86 Workshop Report. In: Schlossman SF, Boumsell L, Gilks W, et al, ed. *Leukocyte Typing V: White Cell Differentiation Antigens*. New York: Oxford University Press; 1995. (Biology)

Jacobsen CN, Aasted B, Broe MK, Petersen JL. Reactivities of 20 anti-human monoclonal antibodies with leucocytes from ten different animal species. *Vet Immunol Immunopathol.* 1993; 39(4):461-466. (Biology)

Schlossman S, Boumsell L, et al, ed. *Leukocyte Typing V*. New York: Oxford University Press; 1995. (Clone-specific)

Wilson AD, Shooshtari M, Finerty S, Watkins P, Morgan AJ. Selection of monoclonal antibodies for the identification of lymphocyte surface antigens in the New World primate *Saguinus oedipus oedipus* (cotton top tamarin). *J Immunol Methods.* 1995; 178(2):195-200. (Biology)