

**Important!**

Includes new buffers!  
Review instructions  
before use.

# ***Pst* I**

Cat. no. 15215-015

Size: 3,000 Units

Store at  $-20^{\circ}\text{C}$

10X Buffer H: 1 mL

10X Loading Buffer: 1 mL

Cleavage Site:

5' -C TGCA↓G-3'

3' -G↑ACGT C-5'

## **Restriction Enzyme Reaction Mix**

*Pst* I 1  $\mu\text{L}$

10X Buffer H 2  $\mu\text{L}$

Substrate DNA  $\leq 1$   $\mu\text{g}$

Sterile water to 20  $\mu\text{L}$

**Incubate at  $37^{\circ}\text{C}$ .**

To use the loading buffer, add 1/10 volume of 10X Loading Buffer to an aliquot of the reaction mix, load on to an agarose gel, and analyze by electrophoresis.

Visit [www.invitrogen.com/restrictionenzymes](http://www.invitrogen.com/restrictionenzymes) for details.

## **Intended Use**

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

**Note:** SDS in the loading buffer may precipitate during storage at room temperature. Warm the buffer to dissolve any SDS precipitate before use.

### Source

*Escherichia coli* ED8654 carrying the plasmid encoding the *Pst* I gene

### Unit Definition

One unit is the amount of enzyme required to completely digest 1  $\mu$ g of  $\lambda$ DNA in 50  $\mu$ L of the reaction mixture in 1 hour at 37°C.

### Relative Activity in Universal Buffers

Buffer	L	M	H	K	T (+BSA)
Relative Activity (%)	<20*	60*	100	80	20*

\*Weak star activity is detected.

### Inactivation by Heat

Enzyme is inactivated by heating at 80°C for 20 minutes.

### Effect of DNA Methylation

Enzyme is not sensitive to dam, dcm, or CG methylation.

### Star Activity

Unrelated site may be cut in the presence of high glycerol concentrations.

## **Buffer Compositions**

### **10X Buffer H**

500 mM Tris-HCl, pH 7.5

100 mM MgCl<sub>2</sub>

10 mM Dithiothreitol

1,000 mM NaCl

### **Storage Buffer**

10 mM Tris-HCl, pH 7.5

100 mM NaCl

10 mM MgCl<sub>2</sub>

0.1 mM EDTA

1 mM DTT

0.15% Triton<sup>®</sup> X-100

0.01% BSA

50% (v/v) glycerol

### **Loading Buffer**

1% SDS

50% glycerol

0.05% Bromophenol blue

## **Product Qualification**

Certificate of Analysis is available from

[www.invitrogen.com/support](http://www.invitrogen.com/support). Search for the certificate of analysis by product lot number, which is printed on the box.

## Related Products

Product	Quantity	Cat. no.
T4 DNA Ligase	250 units	15224-041
T4 Polynucleotide Kinase	1,000 U	18004-028
UltraPure™ Low Melting Point Agarose	100 g	16520-100
TrackIt™ 1 Kb Plus DNA Ladder	100 applications	10488-085
SYBR® Safe DNA Gel Stain (10,000X)	400 µL	S33102
E-Gel® EX Starter Kit, 1%	1 kit	G6511ST

## Technical Support

For technical support, email [tech\\_support@invitrogen.com](mailto:tech_support@invitrogen.com) or visit [www.invitrogen.com](http://www.invitrogen.com)

©2010 Life Technologies Corporation. All rights reserved.  
Carlsbad, CA [www.invitrogen.com](http://www.invitrogen.com) Made in China

The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.  
Triton® X-100 is a registered trademark of Union Carbide Chemicals & Plastics Technology Corp.