

Technical Data Sheet

Human Naive/Memory T Cell Panel

Product Information

Material Number: 561438
Size: 50 Tests
Reactivity: QC Testing: Human

Component:

Description:

Size:

Vol. per Test:

Clone Name:

Storage Buffer:

51-9007097

Alexa Fluor® 647 Mouse Anti-Human CD197 (CCR7)

60 Tests (1 ea)

5 µl

150503

Aqueous buffered solution containing BSA, protein stabilizer, and ≤0.09% sodium azide.

Component:

Description:

Size:

Vol. per Test:

Clone Name:

Storage Buffer:

51-9007098

APC-H7 Mouse Anti-Human CD3 (CD3-epsilon; CD3E; Leu4; T-cell surface antigen T3/Leu-4 epsilon chain; T3E)

60 Tests (1 ea)

5 µl

SK7 (also known as Leu-4)

Aqueous buffered solution containing BSA, protein stabilizer, and ≤0.09% sodium azide.

Component:

Description:

Size:

Vol. per Test:

Clone Name:

Storage Buffer:

51-9007099

PerCP-Cy™5.5 Mouse Anti-Human CD4 (L3T4; Leu3a; T-cell surface antigen T4/Leu-3 ; W3/25 ; CD4 antigen (p55))

60 Tests (1 ea)

5 µl

SK3 (also known as Leu3a)

Aqueous buffered solution containing BSA and ≤0.09% sodium azide.

Component:

Description:

Size:

Vol. per Test:

Clone Name:

Storage Buffer:

51-9007100

FITC Mouse Anti-Human CD45RA

60 Tests (1 ea)

5 µl

HI100

Aqueous buffered solution containing BSA and ≤0.09% sodium azide.

Description

Note: The panel is 50 test size. Extra material for each component for an additional 10 tests is provided for instrument set up purposes.

Multicolor immunofluorescent staining followed by flow cytometric analysis and sorting has led to the phenotypic and functional characterization of multiple peripheral T cell subsets. Three major T cell subsets have been well characterized, i.e., naïve, memory and effector T cells. Naïve T cells have a relatively homogenous cell surface phenotype. They express high levels of CD45RA and the CD197 (CCR7 chemokine receptor). Effector T cells coexpress variable to low levels of CD45RA and no CD197 whereas memory T cells coexpress variable

BD Biosciences

bdbiosciences.com

United States	Canada	Europe	Japan	Asia Pacific	Latin America/Caribbean
877.232.8995	866.979.9408	32.2.400.98.95	0120.8555.90	65.6861.0633	55.11.5185.9995

For country contact information, visit bdbiosciences.com/contact

Conditions: The information disclosed herein is not to be construed as a recommendation to use the above product in violation of any patents. BD Biosciences will not be held responsible for patent infringement or other violations that may occur with the use of our products. Purchase does not include or carry any right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of Becton, Dickinson and Company is strictly prohibited.


For Research Use Only. Not for use in diagnostic or therapeutic procedures. Not for resale.

Unless otherwise noted, BD, BD Logo and all other trademarks are property of Becton, Dickinson and Company. © 2015 BD

561438 Rev. 3



to low levels of CD45RA and high levels of CD197. The Human Naive/Memory T Cell Panel contains fluorescent antibodies (each optimized at 5 µl per test) that are specific for the cell surface antigens: CD45RA, CD197, CD4 and CD3. The channel for detecting Phycoerythrin or Phycoerythrin protein-tandem conjugated proteins has been left open for inclusion of other markers. The panel was designed to standardize the multicolor staining and flow cytometric characterization of the three major CD4+ T cell subsets that arise as a consequence of development or clonal expansion and differentiation driven by antigenic stimulation (eg, in response to allergens, infectious disease or vaccination).

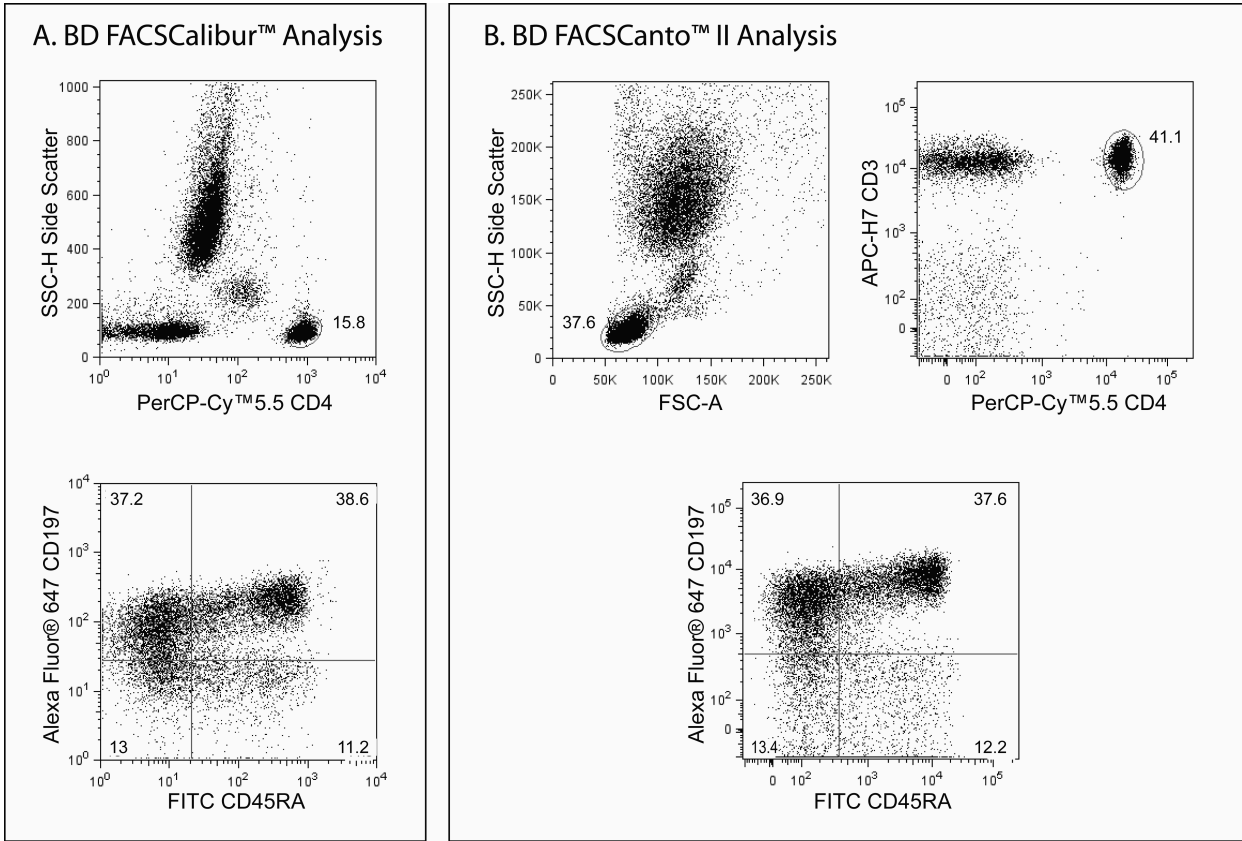

Multicolor staining and flow cytometric analysis of peripheral blood naïve and memory CD4+ T-cells using BD FACSCalibur™ and BD FACSCanto™ Cytometry Systems. Whole blood was stained with a combination of PerCP-Cy™5.5 Mouse anti-Human CD4, FITC Mouse anti-Human CD45RA, and Alexa Fluor® 647 Mouse anti-Human CD197 (CCR7) antibodies for analysis with the BD FACSCalibur™ Cytometer System, or with the addition of APC-H7 Mouse anti-Human CD3 antibody, for analysis with the BD FACSCanto™ Cytometer System. The fluorescent antibodies provided in this kit were each used at 5 µl per test. Red blood cells were then lysed using BD Pharm Lyse™ Lysing Buffer (Cat. No. 555899) and the leukocytes were analyzed. Bivariate dot plots showing the correlated expression patterns of CD45RA versus CD197 (Panels A and B, Lower Plots) were derived from gated events with the light-scattering characteristics of viable CD4+ T lymphocytes (Panel A, Upper Plot, FACSCalibur) or CD4+CD3+ T lymphocytes (Panel B, Upper Plots, FACSCanto II) that were generated by flow cytometric analysis. Quadrants for the dot plots were derived using fluorescence-minus-one (FMO) controls.

Preparation and Storage

Store undiluted at 4°C and protected from prolonged exposure to light. Do not freeze.

The monoclonal antibody was purified from tissue culture supernatant or ascites by affinity chromatography.

The antibody was conjugated to Alexa Fluor® 647 under optimum conditions, and unreacted Alexa Fluor® 647 was removed.

The antibody was conjugated with APC-H7 under optimum conditions, and unconjugated antibody and APC-H7 were removed.

The antibody was conjugated with PerCP-Cy5.5 under optimum conditions, and unconjugated antibody and free PerCP-Cy5.5 were removed. Storage of PerCP-Cy5.5 conjugates in unoptimized diluent is not recommended and may result in loss of signal intensity.

The antibody was conjugated with FITC under optimum conditions, and unreacted FITC was removed.

Application Notes

Application

Flow cytometry	Routinely Tested
----------------	------------------

Suggested Companion Products

Catalog Number	Name	Size	Clone
555899	Lysing Buffer	100 mL	(none)
554656	Stain Buffer (FBS)	500 mL	(none)

BD Biosciences
bdbiosciences.com

United States 877.232.8995	Canada 866.979.9408	Europe 32.2.400.98.95	Japan 0120.8555.90	Asia Pacific 65.6861.0633	Latin America/Caribbean 55.11.5185.9995
--------------------------------------	-------------------------------	---------------------------------	------------------------------	-------------------------------------	---

For country contact information, visit bdbiosciences.com/contact

Conditions: The information disclosed herein is not to be construed as a recommendation to use the above product in violation of any patents. BD Biosciences will not be held responsible for patent infringement or other violations that may occur with the use of our products. Purchase does not include or carry any right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of Becton, Dickinson and Company is strictly prohibited.

For Research Use Only. Not for use in diagnostic or therapeutic procedures. Not for resale.

Unless otherwise noted, BD, BD Logo and all other trademarks are property of Becton, Dickinson and Company. © 2015 BD



Product Notices

1. This reagent has been pre-diluted for use at the recommended Volume per Test. We typically use 1×10^6 cells in a 100- μ l experimental sample (a test).
2. Please refer to www.bdbiosciences.com/pharming/en/protocols for technical protocols.
3. The Alexa Fluor®, Pacific Blue™, and Cascade Blue® dye antibody conjugates in this product are sold under license from Molecular Probes, Inc. for research use only, excluding use in combination with microarrays, or as analyte specific reagents. The Alexa Fluor® dyes (except for Alexa Fluor® 430), Pacific Blue™ dye, and Cascade Blue® dye are covered by pending and issued patents.
4. Alexa Fluor® is a registered trademark of Molecular Probes, Inc., Eugene, OR.
5. Alexa Fluor® 647 fluorochrome emission is collected at the same instrument settings as for allophycocyanin (APC).
6. Caution: Sodium azide yields highly toxic hydrazoic acid under acidic conditions. Dilute azide compounds in running water before discarding to avoid accumulation of potentially explosive deposits in plumbing.
7. For fluorochrome spectra and suitable instrument settings, please refer to our Multicolor Flow Cytometry web page at www.bdbiosciences.com/colors.
8. This APC-conjugated reagent can be used in any flow cytometer equipped with a dye, HeNe, or red diode laser.
9. Please observe the following precautions: Absorption of visible light can significantly alter the energy transfer occurring in any tandem fluorochrome conjugate; therefore, we recommend that special precautions be taken (such as wrapping vials, tubes, or racks in aluminum foil) to prevent exposure of conjugated reagents, including cells stained with those reagents, to room illumination.
10. PerCP-Cy5.5-labelled antibodies can be used with FITC- and R-PE-labelled reagents in single-laser flow cytometers with no significant spectral overlap of PerCP-Cy5.5, FITC, and R-PE fluorescence.
11. PerCP-Cy5.5 is optimized for use with a single argon ion laser emitting 488-nm light. Because of the broad absorption spectrum of the tandem fluorochrome, extra care must be taken when using dual-laser cytometers, which may directly excite both PerCP and Cy5.5™. We recommend the use of cross-beam compensation during data acquisition or software compensation during data analysis.
12. Cy is a trademark of Amersham Biosciences Limited.
13. BD APC-H7 is a tandem conjugate and an analog of APC-Cy7 with the same spectral properties. It has decreased intensity but it is engineered for greater stability and less spillover in the APC channel and consequently offers better performance than APC-Cy7. It has an absorption maximum of approximately 650 nm. When excited by light from a red laser, the APC fluorochrome can transfer energy to the cyanine dye, which then emits at a longer wavelength. The resulting fluorescent emission maximum is approximately 767 nm. BD recommends that a 750-nm longpass filter be used along with a red-sensitive detector such as the Hamamatsu R3896 PMT. As with APC-Cy7 special filters are required when using APC-H7 in conjunction with APC.
Note: Although our APC-H7 products demonstrate higher lot-to-lot consistency than other APC tandem conjugate products, and every effort is made to minimize the lot-to-lot variation in residual emission from APC, it is strongly recommended that every lot be tested for differences in the amount of compensation required and that individual compensation controls are run for each APC-H7 conjugate.
Note: Cy is a trademark of Amersham Biosciences Limited.
14. Although BD APC-H7 is engineered to minimize spillover to the APC channel and is more stable and less affected by light, temperature, and formaldehyde-based fixatives, compared to other APC-cyanine tandem dyes, it is still good practice to minimize as much as possible, any light, temperature and fixative exposure when working with all fluorescent conjugates.
15. Source of all serum proteins is from USDA inspected abattoirs located in the United States.

References

- Appay V, van Lier RA, Sallusto F, Roederer M. Phenotype and function of human T lymphocyte subsets: consensus and issues. *Cytometry A*. 2008; 73(11):975-983. (Biology)
- Kallies A. Distinct regulation of effector and memory T-cell differentiation. *Immunol Cell Biol*. 2008; 86(4):325-332. (Biology)
- Sallusto F, Geginat J, Lanzavecchia A. Central memory and effector memory T cell subsets: function, generation, and maintenance. *Annu Rev Immunol*. 2004; 22:745-763. (Biology)
- Saule P, Trauet J, Dutriez V, Lekeux V, Dessaint JP, Labalette M. Accumulation of memory T cells from childhood to old age: central and effector memory cells in CD4(+) versus effector memory and terminally differentiated memory cells in CD8(+) compartment. 2006; 127(3):274-281. (Biology)
- Seder RA, Darrah PA, Roederer M. T-cell quality in memory and protection: implications for vaccine design. *Nat Rev Immunol*. 2008; 8(4):247-258. (Biology)

BD Biosciences

bdbiosciences.com

United States	Canada	Europe	Japan	Asia Pacific	Latin America/Caribbean
877.232.8995	866.979.9408	32.2.400.98.95	0120.8555.90	65.6861.0633	55.11.5185.9995

For country contact information, visit bdbiosciences.com/contact

Conditions: The information disclosed herein is not to be construed as a recommendation to use the above product in violation of any patents. BD Biosciences will not be held responsible for patent infringement or other violations that may occur with the use of our products. Purchase does not include or carry any right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of Becton, Dickinson and Company is strictly prohibited.

For Research Use Only. Not for use in diagnostic or therapeutic procedures. Not for resale.

Unless otherwise noted, BD, BD Logo and all other trademarks are property of Becton, Dickinson and Company. © 2015 BD

561438 Rev. 3

