

# LIVE/DEAD® Cell Imaging Kit quick start guide

life

kit components (R37601)	amount
Live Green (A) 1 X 10	1mL of 1µM solution
Dead Red (B) 1 X 10	1mM lyophilized stock



1. Thaw Vials



2. Transfer Live Green(A) into Dead Red(B)



3. Mix to create 2X stock



4. Add equal volume 2X stock to cells  
within 2 hours



5. Incubate 15 min at 20-25°C



6. Image cells

**Important Licensing Information** - These products may be covered by one or more Limited Use Label Licenses. By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses.

For research use only. Not for human or animal therapeutic or diagnostic use.

© 2011 Life Technologies Corporation. **All rights reserved.** The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners. R37601\_QuickReferenceCard