

APO-BrdU™ TUNEL Assay Kit

Table 1. Contents and storage information.

Material	Amount	Storage*	Stability
A35125 APO-BrdU™ TUNEL Assay Kit (A23210) ≤-20°C Components			
Positive control cells (Component A, brown cap), fixed human lymphoma cell line†	5 mL	<ul style="list-style-type: none"> • -20°C • Protect from light 	When stored as directed the kit components are stable for at least 3 months.
Negative control cells (Component B, white cap), fixed human lymphoma cell line†	5 mL		
Terminal deoxynucleotidyl transferase (Component C, yellow cap)	45 µL		
5-Bromo-2'-deoxyuridine 5'-triphosphate (BrdUTP) (Component D, violet cap)	480 µL		
A35126 APO-BrdU™ TUNEL Assay Kit (A23210) 2-6°C Components			
Anti-BrdU mouse monoclonal antibody PRB-1, Alexa Fluor 488 conjugate (Component E, orange cap)‡	350 µL	<ul style="list-style-type: none"> • 2-6°C • Protect from light • DO NOT FREEZE 	When stored as directed the kit components are stable for at least 3 months.
Propidium iodide/RNase A staining buffer (Component F, amber bottle)	30 mL		
Reaction buffer (Component G, green cap)§	0.6 mL		
Wash buffer (Component H, blue cap)‡	120 mL		
Rinse buffer (Component I, red cap)‡	120 mL		

*Each module can be stored under the conditions listed. For optimal storage conditions of individual components, refer to the labels on the vials. †The positive and negative control cells are suspensions of $1-2 \times 10^6$ cells/mL in 70% ethanol. ‡The antibody, wash buffer, and rinse buffer each contain 0.05% sodium azide as a preservative. Handle these materials with care. §The reaction buffer contains $\geq 0.2\%$ cacodylic acid (dimethylarsenic) as a buffer. Handle this material with care.

Number of assays: 60 assays of 1 mL samples containing $1-2 \times 10^6$ cells/mL

Approximate fluorescence excitation and emission maxima: 495/519 nm for the Alexa Fluor® 488 labeled monoclonal antibody.

Introduction

A landmark of cellular self-destruction by apoptosis is the activation of nucleases that eventually degrade the nuclear DNA into fragments of approximately 200 base pairs in length.¹ Detection of these DNA fragments is relatively straightforward, making this assay among the most reliable methods for identifying apoptotic cells.² The APO-BrdU™ TUNEL Assay Kit, offered in collaboration with Phoenix Flow Systems, Inc., detects the DNA fragmentation of apoptotic cells by exploiting the fact that the DNA breaks expose a large number

of 3'-hydroxyl ends. These hydroxyl groups can then serve as starting points for terminal deoxynucleotidyl transferase (TdT), which adds deoxyribonucleotides in a template-independent fashion. Addition of the deoxythymidine analog 5-bromo-2'-deoxyuridine 5'-triphosphate (BrdUTP) to the TdT reaction serves to label the break sites. Once incorporated into the DNA, BrdU can be detected by an anti-BrdU antibody using standard immunohistochemical techniques. This method of labeling DNA breaks is referred to as Terminal Deoxynucleotide Transferase dUTP Nick End Labeling, or TUNEL.

The APO-BrdU™ TUNEL Assay Kit provides the materials necessary to label DNA strand breaks for the detection of apoptotic cells, as well as fixed samples of positive and negative control cells for assessing assay performance. Complete protocols for flow cytometry applications are also included, although the kit may be adapted for use with fluorescence microscopy as well. Final detection of BrdU incorporation at DNA break sites is achieved through an Alexa Fluor® 488 dye-labeled anti-BrdU antibody. The Alexa Fluor® 488 dye, which has excitation and emission maxima similar to those of fluorescein, produces fluorescent conjugates that are brighter and more photostable than commonly used fluorescein conjugates. The kit also includes propidium iodide to determine the total cellular DNA content.

Experimental Protocol

The following procedure was designed for use in flow cytometry applications, but is adaptable to fluorescence microscopy [note A]. This protocol requires that cells be in suspension; adherent cell lines may be used but must be trypsinized after induction of apoptosis [note B].

Cell Preparation and Fixation

Cell fixation using paraformaldehyde is a required step in the APO-BrdU™ assay [note C]. The following cell fixation procedure is a suggested method. Variables such as cell origin and growth conditions can affect the results, so the fixation conditions described below should be considered as guidelines. The positive and negative control cells provided in the APO-BrdU™ Kit are already fixed.

- 1.1 Induce apoptosis in cells using the desired method. Although negative control cells are provided with the kit, it may be desirable to prepare a negative control sample using the cell line of interest by incubating cells in the absence of inducing agent.
- 1.2 Suspend $1-2 \times 10^6$ cells in 0.5 mL of phosphate-buffered saline (PBS).
- 1.3 Add the cell suspension into 5 mL of 1% (w/v) paraformaldehyde in PBS and place on ice for 15 minutes.
- 1.4 Centrifuge the cells for 5 minutes at $300 \times g$ and discard the supernatant.
- 1.5 Wash the cells in 5 mL of PBS then pellet the cells by centrifugation. Repeat.
- 1.6 Resuspend the cells in 0.5 ml of PBS.
- 1.7 Add the cells to 5 mL of ice-cold 70% (v/v) ethanol. Let the cells stand for a minimum of 30 minutes on ice or in a -20°C freezer. In some biological systems, storage of the cells at -20°C in 70% (v/v) ethanol for at least 12–18 hours prior to performing the TUNEL assay yields the best results. Cells can be stored at -20°C for several days before use.

Detection of Apoptotic Cells

The following protocol describes the method for measuring apoptosis in the positive and negative control cells that are provided in the APO-BrdU™ TUNEL Assay Kit. The same procedure should be employed for measuring apoptosis in the experimental samples.

- 2.1 Resuspend the positive (brown cap) and negative (white cap) control cells by swirling the vials. Remove 1 mL aliquots of the control cell suspensions (approximately 1×10^6 cells/mL) and place in 12×75 mm flow cytometry centrifuge tubes [note D]. Centrifuge ($300 \times g$) the control cell suspensions for 5 minutes and remove the 70% (v/v) ethanol by aspiration, being careful to not disturb the cell pellet.
- 2.2 Resuspend the control cells of each tube with 1 mL of wash buffer (blue cap). Centrifuge for 5 minutes at $300 \times g$ and remove the supernatants by aspiration. Repeat.
- 2.3 Prepare a DNA-labeling solution; a total volume of 50 μ L is required for each sample. Mix 10 μ L of reaction buffer (green cap), 0.75 μ L of TdT enzyme (yellow cap), 8.0 μ L of BrdUTP (violet cap) and 31.25 μ L of dH₂O. For additional samples the volumes may be scaled up accordingly, but mix only enough DNA-labeling solution to complete the number of assays prepared per session. The DNA-labeling solution is active for approximately 24 hours.
- 2.4 Resuspend the control cell pellets of each tube in 50 μ L of the DNA-labeling solution (prepared in step 2.3).
- 2.5 Incubate the cells in the DNA-labeling solution for 60 minutes at 37°C in a temperature controlled bath. Shake the samples every 15 minutes to keep the cells in suspension. For samples other than the control cells provided in the kit, incubation times at 37°C may need to be adjusted to longer or shorter periods depending on the characteristics of the experimental samples. The DNA-labeling reaction for the control cells can also be carried out at 22–24°C overnight.
- 2.6 At the end of the incubation time add 1.0 mL of rinse buffer (red cap) to each tube and centrifuge at $300 \times g$ for 5 minutes. Remove the supernatants by aspiration.
- 2.7 Repeat the cell rinsing (as in step 2.6) with 1.0 mL of rinse buffer (red cap). Centrifuge the samples at $300 \times g$ and remove the supernatants by aspiration.
- 2.8 Prepare 100 μ L of antibody staining solution for each sample by mixing 5.0 μ L of the Alexa Fluor® 488 dye-labeled anti-BrdU antibody (orange cap) with 95 μ L of rinse buffer (red cap). Prepare only enough antibody staining solution to complete the number of assays prepared per session.
- 2.9 Resuspend the cell pellets in 100 μ L of the antibody solution prepared in step 2.8. Incubate the cells in this solution for 30 minutes at room temperature. Protect the samples from light during the incubation.
- 2.10 Add 0.5 mL of the propidium iodide/RNase A staining buffer (amber bottle) to each sample [note E]. Incubate the cells for an additional 30 minutes at room temperature. Protect the samples from light during the incubation.
- 2.11 Analyze the samples by flow cytometry [notes F and G]. It is recommended that the samples be analyzed within 3 hours of completing the staining procedure.

Product Notes

[A] For microscopy applications, it is recommended that the cells be deposited onto slides after the antibody staining step, but prior to the propidium iodide/RNase treatment. Cells that have undergone apoptosis should fluoresce brightly when viewed with filter sets appropriate for fluorescein.

[B] For adherent cell lines, detached cells present in the supernatant have a higher probability of being apoptotic than do cells that have remained adherent. Detached cells should be collected prior to trypsinization of the adherent cell layer.

[C] Cell fixation is an important step in analyzing apoptotic samples. Unfixed cells may lose smaller DNA fragments, leading to lower signals. Alternate fixation methods may be necessary to fully exploit some cell systems.

[D] To minimize cell loss, use a single 12 × 75 mm test tube for each assay. Use of sequential tubes can result in significant cell loss during the assay due to cells sticking to the sides of the tube.

[E] If the DNA cell cycle information is not required, it is not necessary to add the PI/RNase A staining buffer to each tube.

[F] If signal intensity is low, the incubation time for the DNA-labeling reaction (step 2.5) may be extended. Labeling times of up to 4 hours at 37°C may be required for some systems.

[G] Occasionally a mirror image population of cells at lower intensity is observed in the flow cytometry dual parameter display. This population arises when some cells become stuck to the side of the test tube and are not fully exposed to the DNA-labeling solution. Make sure all cells are properly suspended at the beginning of the labeling reaction.

References

1. Am J Pathol 136, 593 (1990); 2. Trends Cell Biol 5, 21 (1995).

Product List

Current prices may be obtained from our website or from our Customer Service Department.

Cat #	Product Name	Unit Size
A23210	APO-BrdU™ TUNEL Assay Kit *with Alexa Fluor® 488 anti-BrdU* *60 assays*.....	1 kit

Contact Information

Molecular Probes, Inc.
29851 Willow Creek Road
Eugene, OR 97402
Phone: (541) 465-8300
Fax: (541) 335-0504

Customer Service:
6:00 am to 4:30 pm (Pacific Time)
Phone: (541) 335-0338
Fax: (541) 335-0305
probesorder@invitrogen.com

Toll-Free Ordering for USA:
Order Phone: (800) 438-2209
Order Fax: (800) 438-0228

Technical Service:
8:00 am to 4:00 pm (Pacific Time)
Phone: (541) 335-0353
Toll-Free (800) 438-2209
Fax: (541) 335-0238
probetech@invitrogen.com

Invitrogen European Headquarters
Invitrogen, Ltd.
3 Fountain Drive
Inchinnan Business Park
Paisley PA4 9RF, UK
Phone: +44 (0) 141 814 6100
Fax: +44 (0) 141 814 6260
Email: euroinfo@invitrogen.com
Technical Services: eurotech@invitrogen.com

Further information on Molecular Probes products, including product bibliographies, is available from your local distributor or directly from Molecular Probes. Customers in Europe, Africa and the Middle East should contact our office in Paisley, United Kingdom. All others should contact our Technical Service Department in Eugene, Oregon.

Molecular Probes products are high-quality reagents and materials intended for research purposes only. These products must be used by, or directly under the supervision of, a technically qualified individual experienced in handling potentially hazardous chemicals. Please read the Material Safety Data Sheet provided for each product; other regulatory considerations may apply.

Limited Use Label License No. 223: Labeling and Detection Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) to not transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of the above patents based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Molecular Probes, Inc., Business Development, 29851 Willow Creek Road, Eugene, OR 97402, Tel: (541) 465-8300. Fax: (541) 335-0354.

Several Molecular Probes products and product applications are covered by U.S. and foreign patents and patents pending. All names containing the designation ® are registered with the U.S. Patent and Trademark Office.

Copyright 2006, Molecular Probes, Inc. All rights reserved. This information is subject to change without notice.