



Qty: 2 ml  
 rec-Protein G-Sepharose® 4B Conjugate  
 Catalog No. 10-1241  
 Lot No.: See product label  
 Exp. Date: See product label

### rec-Protein G-Sepharose® 4B Conjugate

#### FORM

This product is supplied as a 50% suspension containing 2 ml rec-Protein G-Sepharose 4B + 2 ml PBS (10 mM phosphate-buffered saline, pH 7.4, containing 0.1% Proclin preservative). This product is prepared by covalently coupling highly purified rec-Protein G (produced in *E. Coli*) to Sepharose® 4B.

#### DESCRIPTION

Protein A and Protein G bind specifically to Fc regions of many mammalian immunoglobulins. Protein A and Protein G conjugates are commonly used as affinity adsorbents to purify immunoglobulins (antibodies) and immunoglobulin subtypes from serum, hybridoma ascites fluids, tissue culture supernatants and other biological fluids. These reagents are also commonly used to capture immune complexes generated in immunoprecipitation experiments. Because Protein A and Protein G vary widely in their abilities to bind individual immunoglobulins, consult the table below before proceeding with your application. It should be noted that the Fc binding characteristics of naturally produced Protein A and Protein G are identical to their recombinant counterparts. However, also note that Invitrogen's recombinant Protein G products are deleted for two albumin binding sites present in the wild type molecule. For certain procedures, it may be advantageous to use a light chain binding reagent (e.g. Invitrogen's Kappalock™ products) instead of Protein A or Protein G.

#### SPECIFICATIONS

- rec-Protein-G/Sepharose® ratio:** ~2.5 mg rec-Protein G per ml Sepharose
- Binding capacity:** ~18 mg Rabbit IgG per ml beads
- rec-Protein G molecular weight:** 22 kDa (migrates as 32-36 by SDS-PAGE)
- Ig Binding Sites/molecule rec-Protein G:** 2 per molecule
- Albumin binding sites/molecule rec-Protein G:** none

**Binding Characteristics of Immunoglobulins to Protein A and Protein G**

Species	Protein A	Protein G	Species	Protein A	Protein G	Species	Protein A	Protein G
Cat IgG	+++	+	Human IgA2	+	-	Mouse IgG2b	+++	+++
Chicken IgG	-	-	Human IgD	-	-	Mouse IgG3	++	+++
Cow IgG	+	+++	Human IgE	++	-	Mouse IgM	-	-
Cow IgG1	+	+++	Human IgG	+++	+++	Rat IgG	+	++
Cow IgG2	+++	+++	Human IgG1	+++	+++	Rat IgG1	-	+
Dog IgG	+++	+	Human IgG2	+++	+++	Rat IgG2a	-	+++
Goat IgG	+	++	Human IgG3	-	+++	Rat IgG2b	-	+
Goat IgG1	+	+++	Human IgG4	+++	+++	Rat IgG2c	++	++
Goat IgG2	+++	+++	Human IgM	+	-	Rabbit IgG	+++	+++
Guinea Pig IgG	+++	+	Monkey IgG	+++	+++	Sheep IgG	+	++
Hamster IgG	+	++	Pig IgG	+++	++	Sheep IgG1	+	++
Horse IgG	+	+++	Mouse IgG	++	++	Sheep IgG2	+++	+++
Human IgA	+	-	Mouse IgG1	+	++			
Human IgA1	+	-	Mouse IgG2a	+++	+++			

+++ , ++ , + , --: strong binding, moderate binding, weak binding, no binding, respectively.

(cont'd)

**For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.**

[www.invitrogen.com](http://www.invitrogen.com)

Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288 • E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)

PI101241

(Rev 06/10) DCC-10-1592

**Important Licensing Information** - These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or our website, [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses. Unless otherwise indicated, these products are for research use only and are not intended for human or animal diagnostic, therapeutic or commercial use.

**PROCEDURES:**

Recommended procedures for immunoglobulin affinity purification and immunoprecipitation using our Protein A and Protein G products are available on our Web site at [www.invitrogen.com/methods](http://www.invitrogen.com/methods). Hard copy is available on request from Invitrogen Technical Services by phone at 800.955.6288 or by email at [tech\\_support@invitrogen.com](mailto:tech_support@invitrogen.com). An excellent source for information on antibodies in general and on most routine laboratory procedures involving antibodies is Antibodies, A Laboratory Manual, Ed Harlow and David Lane, Eds., Cold Spring Harbor Laboratory (1999).

**STORAGE:**

Store at 2-8°C. **Do Not Freeze.** Freezing this product will damage or destroy its performance.

**SELECTED INVITROGEN PRODUCT REFERENCES***Immunoprecipitation*

1. Dikic I, et al. *Nature* 383(6600):547-550, 1996.
2. Streuli CH, et al. *Cell* 69:927-938, 1995.
3. Goi K, et al. *Cancer Research* 57(10):1895-1902, 1997.
4. Elliott JG, et al. *J Biol Chem* 272(21):13849-13855, 1997.
5. Inohara H, et al. *Cancer Research* 56:4530-4534, 1996.
6. McConnell KR, et al. *J Immunol* 158:2083-2089, 1997.

*Affinity Purification*

7. Sakuai T, et al. *J Cell Biol* 136:907-918, 1997.
8. Lasky L, et al. *Cell* 69:927-938, 1992.

**RELATED PRODUCTS****Product**

Product	Cat. No.
Protein A-Sepharose® (5 ml)	10-1041
Protein A-Sepharose® (25 ml)	10-1142
Rec-Protein A-Sepharose® (25 ml)	10-1142
MonoAb rec-Protein A Column Kit	10-1160
Zysorbin™ (fixed, killed Protein A-bearing <i>Staph. aureus</i> for IP, 25 ml)	10-1051-1
Zysorbin™ (fixed, killed Protein A-bearing <i>Staph. aureus</i> for IP, 100 ml)	10-1051-3
Rec-Protein G-Sepharose® (5 ml)	10-1242
Zysorbin-G (fixed, killed Protein G-bearing Streptococci for IP, 10 ml)	10-1352
Kappalock™-Sepharose (binds to mammalian kappa light chains, 2 ml)	10-1841

Zymed® is a registered trademark of Invitrogen. Sepharose® is a registered trademark of Amersham Biosciences Limited. Zysorbin™ is a Zymed trademark. Kappalock™ is a trademark of Aston Inc.

**Explanation of symbols**

Symbol	Description	Symbol	Description
	Catalogue Number		Batch code
	Research Use Only		In vitro diagnostic medical device
	Use by		Temperature limitation
	Manufacturer		European Community authorised representative
	Without, does not contain		With, contains
	Protect from light		Consult accompanying documents
	Directs the user to consult instructions for use (IFU), accompanying the product.		

**For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.**

[www.invitrogen.com](http://www.invitrogen.com)

Invitrogen Corporation • 542 Flynn Rd • Camarillo • CA 93012 • Tel: 800.955.6288 • E-mail: [techsupport@invitrogen.com](mailto:techsupport@invitrogen.com)

PI101241

(Rev 06/10) DCC-10-1592

**Important Licensing Information** - These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or our website, [www.invitrogen.com](http://www.invitrogen.com)). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses. Unless otherwise indicated, these products are for research use only and are not intended for human or animal diagnostic, therapeutic or commercial use.