

## Secondary (2°) Antibody Solutions Alk-Phos. Conjugates

Cat. nos. WP20006, WP20007

Store at 4°C

### Description

Secondary (2°) antibody solutions are pre-diluted, ready-to-use formulations containing polyclonal antibodies conjugated to alkaline phosphatase (AP) in a proprietary stabilizing reagent. The 2° antibody conjugates are specially purified to reduce cross reactivity with bovine antigens. Available 2° antibody solutions include goat anti-rabbit and goat anti-mouse polyclonal alkaline phosphatase conjugates. Secondary antibody solutions can be used for either chemiluminescent or chromogenic detection of antigens bound with a primary antibody of the appropriate host species.

### Products

Cat. no.	Description	Volume
WP20006	2° Antibody Solution Alk-Phos. Conjugated (Anti-Mouse)	2 × 100 ml
WP20007	2° Antibody Solution Alk-Phos. Conjugated (Anti-Rabbit)	2 × 100 ml

### Storage Conditions

Store all reagents at 4°C. Do not freeze 2° antibody solutions.

### Safety Considerations

There are no hazardous chemicals in these conjugate solutions. However, the formulations do contain dilute solutions of irritants. Wear gloves, safety glasses, and a lab coat when using these reagents.

### Method

An example of a standard western blot or western dot blot detection procedure for nitrocellulose (NC) or polyvinylidene fluoride (PVDF) membranes is described below. The protocol can be modified based on initial results.

### Materials Required

- Blotted membrane with antigen of interest
- Purified water — autoclaved or ultrafiltered to remove alkaline phosphatase activity
- Clean containers for preparing solutions and incubating the membrane
- Clean forceps for manipulating blotted membrane
- Orbital shaker capable of rotating at 1 revolution/second
- Blocking buffer such as WesternBreeze® Blocker/Diluent (see page 4) or 5% non-fat dry milk in Tris Buffered Saline with Tween (0.05 M Tris-HCl, 0.15 M NaCl, 0.1% Tween 20, pH 7.5)
- Wash buffer such as WesternBreeze® Wash Solution (see page 4)
- High affinity antigen-specific primary antibody diluted in blocking buffer at manufacturer recommended concentrations
- 2° Antibody Solution Alk-Phos. Conjugate specific to the primary antibody host animal
- Chromogenic or chemiluminescent AP substrate (see page 4)

### General Guidelines

To obtain the best results with 2° antibody solutions:

- Use enough solution to keep the membrane completely covered at all times.
- Use a single, clean dish for each blot. Containers should be large enough to accommodate the membrane and allow it to be fully immersed by solutions.
- Avoid touching the surface of the membrane. Wear clean gloves and handle the blot only with clean forceps.
- Use pure water, free from alkaline phosphatase activity for making solutions and washes. Fresh ultra-filtered water is preferred. Autoclave or ultra-filter stored water to remove alkaline phosphatase activity.
- Avoid cross-contamination of system solutions with the alkaline phosphatase substrate solution.
- Perform all washing, blocking, and incubation steps on a rotary orbital shaker rotating at 1 revolution/second.
- Work quickly when changing solutions to ensure membranes remain wet. PVDF membranes dry quickly and must be re-wet with methanol and rinsed with water before proceeding if they become dry (see **Preparing Stored Membranes**).

### General Guidelines, Continued

- Add solutions to the trays slowly, at the membrane edge, to avoid bubble formation under the membrane. Decant from the same corner of the dish to ensure complete removal of previous solutions.

### Suggestions for Antibody Usage

Antibody concentrations for chemiluminescent detection are lower than those used for chromogenic detection. For best result empirically determine the optimal concentration of antibody needed by dot blot. Antibody solutions that are too dilute result in weak or no signal, whereas overly concentrated solutions cause high background or non-specific binding. Dilute primary antibodies in blocking buffer (see **Materials Required**). **Do not** dilute 2° antibody solutions.

### Preparing Stored Membranes

Membranes can be stored after transfer for subsequent immunodetection after being washed to remove gel and transfer buffer components. When starting with washed and dried membranes, perform the following steps:

Membrane	Procedure
NC	Wash the NC membrane with water for 5 minutes twice, then proceed directly to step 2 of the <b>Western Blot Procedure</b> .
PVDF	Re-wet the PVDF membrane in methanol and then wash with water for 5 minutes twice before proceeding to step 2 of the <b>Western Blot Procedure</b> .

When performing native-PAGE western blots, a drying step is recommended to improve protein binding to the membrane. To re-wet the membrane, follow the appropriate instructions from the table above.

### Western Blot Procedure

1. After transfer, wash the membrane with water for 5 minutes twice, to remove any gel and transfer buffer components.
2. Place the membrane in blocking buffer at room temperature for 10–60 minutes with shaking (1 revolution/sec), or if desired, overnight at 4°C without shaking. Decant blocking buffer.
3. Add diluted primary antibody solution. Incubate for 30–60 minutes with shaking. Decant solution.
4. Add an excess volume of wash buffer. Wash for 1 minute with shaking. Decant solution.
5. Add a volume of fresh wash buffer and wash with shaking for 5 minutes, then decant. Repeat this step up to 3 times.
6. Add pre-diluted, ready-to-use 2° Antibody Solution. Incubate for 30 minutes. Decant solution.
7. Add an excess volume of wash buffer. Wash for 1 minute with shaking. Decant solution.
8. Add another volume of wash buffer and wash with shaking for 5 minutes, then decant. Repeat this step up to 3 times.  
**Note: Do not use phosphate buffer to wash.** Inorganic phosphate is a powerful inhibitor of AP.
9. Rinse the blot briefly with water for 2 minutes, then decant. Proceed immediately to detection.

### Chemiluminescent Detection

1. For chemiluminescent detection, prepare substrate according to manufacturer's instructions.
2. Lay the blot on a clean piece of plastic sheet (such as clear copier transparency material), antigen side up. **Do not** allow the membrane to dry.
3. Pipet the chemiluminescent substrate working solution onto the blot surface, taking care to cover the entire blot.
4. Allow the reaction to develop for the amount of time recommended by the manufacturer.
5. Absorb excess substrate by gently blotting with clean absorbent material (such as filter paper or tissue) or by tilting the plastic sheet so excess substrate drains off onto the absorbent material, but **do not** let the membrane dry.
6. Overlay another clean piece of plastic on top of the blot to prevent drying during imaging. Gently lowering the top plastic from one side of the blot to the other will prevent bubble formation on top of the blot. If excess substrate squeezes out from between the plastic sheets take care to blot it away. If X-ray film imaging is used, no liquid must contact the film. Plastic wrap can be used as an alternative to plastic sheets.
7. Place the protected blot in a film cassette or imaging system and expose for 1 second to several minutes depending on the signal produced and image desired.

### Chromogenic Detection

1. For chromogenic detection, add enough chromogenic substrate to cover the blot. Incubate with shaking until the desired band intensity is achieved. Decant solution.
2. Stop reaction by rinsing briefly with water and decanting the wash. Rinse with water for 2 minutes. Repeat 2 minute water rinse twice.
3. Dry the membrane on clean filter paper to open air or under an infrared lamp.
4. Record an image of the blot. Store membrane protected from light to prevent band fading. Bands remain visible for years when protected from light.

## Troubleshooting

Review the information below to troubleshoot your experiments. For additional troubleshooting, refer to the manual for the WesternBreeze® Chromogenic Western Blot Immunodetection or WesternBreeze® Chemiluminescent Western Blot Immunodetection Kit available at [www.invitrogen.com](http://www.invitrogen.com).

<b>Problem</b>	<b>Possible cause</b>	<b>Possible Solution</b>
<b>Weak or No Signal</b>	Poor or incomplete transfer	Make sure transfer apparatus and membrane sandwiches are assembled correctly. Use appropriate transfer times. Repeat blot. After blotting, stain membrane to measure transfer efficiency.
	Protein of interest ran off the gel	Use positive control and/or molecular weight marker to match gel separation range to size of protein being blotted. After blotting, stain membrane to measure transfer efficiency.
	Membrane not completely wetted	Follow instructions for pre-wetting the membrane.
	Inactive or overly dilute primary antibody	Determine antibody activity by performing a dot blot. Increase antibody concentration as necessary.
	Sample improperly prepared or antigenicity compromised	SDS and reducing agents may interfere with some antibody/antigen affinities.
	Sample too dilute	Load a larger amount of protein onto the gel.
	Poor retention of proteins or protein weakly bound to membrane	Ensure that transfer buffer contains 10–20% methanol. Use membranes with appropriate binding capacity.
	Insufficient exposure time (chemiluminescent detection)	Re-expose film for a longer period of time.
	Insufficient substrate incubation (chromogenic detection)	Remove blot from substrate when signal-to-noise ratio is acceptable.
	Blots are too old	Protein may have degraded over time. Use freshly prepared blots.
<b>High Background</b>	Film overexposed or became wet during exposure (chemiluminescent detection)	Decrease exposure time or allow signal to further decay. Prevent leakage of solutions by encasing membrane in transparency film and blotting excess substrate from edges before exposure.
	Blot is overdeveloped (chromogenic detection)	Follow recommended developing time or remove blot from substrate when signal-to-noise ratio is acceptable.
	Blocking time or washing time is too short	Perform each step for the specified amount of time.
	Primary antibody concentration too high	Determine optimal antibody concentration by performing a dot blot. Decrease antibody concentration as necessary.
	Membrane not completely wetted	Follow instructions for pre-wetting the membrane. The incubation dish must be small enough to allow thorough coverage of membrane to prevent drying. Shake or agitate during each step.
	Membrane, solutions, or incubation tray is contaminated	Use clean glassware and purified water to prepare solutions. Replace or clean the tray thoroughly with a glassware-cleaning detergent. Rinse thoroughly with purified water. Wear clean gloves at all times. Use forceps when handling membranes.
<b>Non-Specific Binding</b>	Insufficient removal of SDS or weakly bound proteins from membrane after blotting	Follow instructions for membrane preparation before immunodetection.
	Short blocking time or long washing time	Make sure that each step is performed for the specified amount of time.
	Affinity of the primary antibody for the protein standards	Check with protein standard manufacturer for homologies with primary antibody.
	Membrane is contaminated by fingerprints or keratin proteins	Wear clean gloves at all times and use forceps when handling membranes. Always handle membranes around the edges.

## Troubleshooting, Continued

<b>"Spotted" Membrane</b>	Membrane blotting pads are dirty or contaminated	Soak pads with detergent and rinse thoroughly with purified water before use. Replace pads when they become worn or discolored.
	Blocking was uneven	The incubation dish must be sufficient to allow thorough coverage of membrane. Shake or agitate during each step.
<b>Large, Scattered Signal</b>	Protein is overloaded	Reduce load or dilute concentration of sample.
	Poor or incomplete transfer	Make sure transfer apparatus and membrane sandwiches are assembled correctly. Use appropriate transfer times. Repeat blot.

## Related Products

Product	Amount	Catalog no.
WesternBreeze® Blocker/Diluent (part A and B)	2 × 80 ml	WB7050
WesternBreeze® Wash Solution (16X)	2 × 100 ml	WB7003
Novex® AP Chromogenic Substrate	250 ml	WP20001
Novex® AP Chemiluminescent Substrate	100 ml	WP20002
Novex® AP Chemiluminescent Substrate Enhancer	5 ml	WP20003
iBlot™ Gel Transfer Device	1 unit	IB1001
iBlot™ Transfer Stack, Regular (Nitrocellulose)	10 sets/box	IB3010-01
iBlot™ Transfer Stack, Mini (Nitrocellulose)	10 sets/box	IB3010-02
iBlot™ Transfer Stack PVDF, Regular	10 sets/box	IB4010-01
iBlot™ Transfer Stack PVDF, Mini	10 sets/box	IB4010-02
Nitrocellulose, 0.2 µm/Filter Paper Sandwiches	20 membranes	LC2000
Nitrocellulose, 0.45 µm/Filter Paper Sandwiches	20 membranes	LC2001
PVDF, 0.2 µm/Filter Paper Sandwiches	20 membranes	LC2002
Invitrolon™ PVDF/Filter Paper Sandwiches	20 membranes	LC2005
Novex® Reversible Membrane Protein Stain Kit	1 kit	IB7710
Sponge Pad for blotting	8 pads	EI9052
Primary antibody	---	visit <a href="http://www.invitrogen.com/antibodies">www.invitrogen.com/antibodies</a>

## Product Qualification

The Certificate of Analysis (CofA) provides detailed quality control information for each product. The CofA is available on our website at [www.invitrogen.com/cofa](http://www.invitrogen.com/cofa), and is searchable by product lot number, which is printed on each box.

## Limited Use Label License No. 5: Invitrogen Technology

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing (a) not to transfer such materials to any third party, and (b) to use such transferred materials and/or information solely for research and not for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. Invitrogen Corporation will not assert a claim against the buyer of infringement of patents owned or controlled by Invitrogen Corporation which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product. If the purchaser is not willing to accept the limitations of this limited use statement, Invitrogen is willing to accept return of the product with a full refund. For information on purchasing a license to this product for purposes other than research, contact Licensing Department, Invitrogen Corporation, 5791 Van Allen Way, Carlsbad, California 92008. Phone (760) 603-7200. Fax (760) 602-6500. Email: [outlicensing@invitrogen.com](mailto:outlicensing@invitrogen.com).

©2008 Invitrogen Corporation. All rights reserved.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.