



ELISA Kit
Catalog #KHR9121

**Insulin
Receptor*
[pY1158]**

www.invitrogen.com

Invitrogen Corporation

542 Flynn Road, Camarillo, CA 93012

Tel: 800-955-6288

E-mail: techsupport@invitrogen.com

*Patent Pending

TABLE OF CONTENTS

Introduction	4
Principle of the Method.....	6
Reagents Provided.....	7
Supplies Required But Not Provided	8
Procedural Notes/Lab Quality Control.....	8
Safety	10
Directions for Washing.....	10
Procedure for Extraction of Proteins from Cells.....	11
Reagent Preparation and Storage.....	13
Reconstitution and Dilution of IR [pY1158] Standard	13
Dilution of IR [pY1158] Standard.....	14
Storage and Final Dilution of Anti-Rabbit IgG HRP (100X).....	15
Dilution of Wash Buffer	16
Assay Method.....	16
Typical Data	19
Limitations of the Procedure.....	20
Performance Characteristics	21
Sensitivity	21
Precision	22
Recovery.....	23
Parallelism	23
Linearity of Dilution	24
Specificity.....	25
References	28

INTRODUCTION

Insulin receptor (IR), a cell surface receptor, binds insulin and mediates its action on target cells. Insulin receptor belongs to the superfamily of the growth factor receptor tyrosine kinases that regulate multiple signaling pathways through activation of a series of phosphorylation cascades. The insulin receptor is a heterotetrameric membrane glycoprotein consisting of disulfide-linked subunits in a β - α - α - β configuration. The α -subunit (135 kDa) is completely extracellular, whereas the β -subunit (95 kDa) possesses a single transmembrane domain with tyrosine kinase activity. Insulin binding to the extracellular domain leads to autophosphorylation of the receptor and activation of the intrinsic tyrosine kinase activity, which allows appropriate substrates to be phosphorylated.

Once activated, the IR initiates a variety of metabolic functions including glucose transport, glycogen synthesis, protein synthesis, translational control and mitogenesis. The IR also plays an important role in neurological function and in hypertension. Important down-stream proteins associated with IR signaling include IRS-1, JAK1, JAK2, STAT1, STAT3, PI3 kinase, GSK-3 β , PKC, SHC, ERK1/2 and many others.

Insulin receptor and insulin-like growth factor 1 receptor (IGF-1R), which are structurally related, share conserved tyrosine residues that are phosphorylated in IR in response to insulin and in IGF-1R in response to IGF-1. Functionally, the insulin receptor regulates metabolism and IGF-1R mediates growth and differentiation. The catalytic loops within the tyrosine kinase domains of the IR/IGF-1R share the same residue sequence with a three-tyrosine motif corresponding to Tyr1158, 1162

and 1163 (for the IR) and Tyr1131, 1135 and 1136 (for the IGF-1R). It is generally believed that autophosphorylation within the activation loop proceeds in a progressive manner initiating at the second tyrosine (1162 or 1135), followed by phosphorylation at the first tyrosine (1158 or 1131), then the last (1163 or 1136), upon which the IR or IGF-1R becomes fully active.

The Invitrogen IR [pY1158] ELISA is designed to detect and quantify the levels of insulin receptor β -subunit that are phosphorylated at residue tyrosine 1158 of insulin receptor. Both natural (heterotetrameric) and recombinant IR react in this assay. This ELISA does not detect IGF-R phosphorylated at tyrosine 1131. This assay allows differentiation of the phosphorylation and activation of IR from that of IGF-1R. Although this ELISA kit is developed using human cells, cross-reactivity with mouse and rat insulin receptor is documented. This assay is intended for detection of IR [pY1158] from lysates of cells. For normalizing the IR content of the samples, an IR (β -subunit) ELISA kit which is independent of phosphorylation status is available from Invitrogen (Cat. #KHR9111).

For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.

READ ENTIRE PROTOCOL BEFORE USE

PRINCIPLE OF THE METHOD

The Invitrogen Insulin Receptor [pY1158] kit is a solid phase sandwich Enzyme Linked-Immuno-Sorbent Assay (ELISA). A monoclonal antibody specific for IR (β -subunit) (regardless of phosphorylation state) has been coated onto the wells of the microtiter strips provided. Samples, including a standard containing IR [pY1158], control specimens and unknowns, are pipetted into these wells. During the first incubation, the IR antigen binds to the immobilized (capture) antibody. After washing, an antibody specific for IR phosphorylated at tyrosine 1158 is added to the wells. During the second incubation, this antibody serves as a detection antibody by binding to the immobilized IR protein captured during the first incubation. After removal of excess detection antibody, a horseradish peroxidase-labeled Anti-Rabbit IgG (Anti-Rabbit IgG HRP) is added. This binds to the detection antibody to complete the four-member sandwich. After a third incubation and washing to remove all the excess Anti-Rabbit IgG HRP, a substrate solution is added, which is acted upon by the bound enzyme to produce color. The intensity of this colored product is directly proportional to the concentration of IR [pY1158] present in the original specimen.

REAGENTS PROVIDED

Note: Store all reagents at 2 to 8°C.

<i>Reagents</i>	<i>96 Test Kit</i>
<i>IR [pY1158] Standard:</i> Contains 0.1% sodium azide. Refer to vial label for quantity and reconstitution volume.	2 vials
<i>Standard Diluent Buffer:</i> Contains 0.1% sodium azide; 25 mL per bottle.	1 bottle
<i>Antibody Coated Wells.</i> 12 x 8 Well Strips.	1 plate
<i>IR [pY1158] Detection Antibody.</i> Contains 0.1% sodium azide; 11 mL per bottle.	1 bottle
<i>Anti-Rabbit IgG HRP (100X).</i> Contains 3.3 mM thymol; 0.125 mL per vial.	1 vial
<i>HRP Diluent.</i> Contains 3.3 mM thymol; 25 mL per bottle.	1 bottle
<i>Wash Buffer Concentrate (25X).</i> 100 mL per bottle.	1 bottle
<i>Stabilized Chromogen, Tetramethylbenzidine (TMB).</i> 25 mL per bottle.	1 bottle
<i>Stop Solution.</i> 25 mL per bottle.	1 bottle
<i>Plate Covers, adhesive strips.</i>	3

Disposal Note: This kit contains materials with small quantities of sodium azide. Sodium azide reacts with lead and copper plumbing to form explosive metal azides. Upon disposal, flush drains with a large volume of water to prevent azide accumulation. Avoid ingestion and contact with eyes, skin and mucous membranes. In case of contact, rinse affected area with plenty of water. Observe all federal, state and local regulations for disposal.

SUPPLIES REQUIRED BUT NOT PROVIDED

1. Microtiter plate reader capable of measurement at or near 450 nm.
2. Calibrated adjustable precision pipettes, preferably with disposable plastic tips. (A manifold multi-channel pipette is desirable for large assays.)
3. Cell extraction buffer (see Recommended Formulation, p. 11).
4. Distilled or deionized water.
5. Plate washer: automated or manual (squirt bottle, manifold dispenser, etc.).
6. Data analysis and graphing software. Graph paper: linear (Cartesian), log-log, or semi-log, as desired.
7. Glass or plastic tubes for diluting and aliquoting standard.
8. Absorbent paper towels.
9. Calibrated beakers and graduated cylinders in various sizes.

PROCEDURAL NOTES/LAB QUALITY CONTROL

1. When not in use, kit components should be refrigerated. All reagents should be warmed to room temperature before use.
2. **Microtiter plates should be allowed to come to room temperature before opening the foil bags.** Once the desired number of strips has been removed, immediately reseal the bag and store at 2 to 8°C to maintain plate integrity.

3. Samples should be frozen if not analyzed shortly after collection. Avoid multiple freeze-thaw cycles of frozen samples. Thaw completely and mix well prior to analysis.
4. If particulate matter is present, centrifuge or filter prior to analysis.
5. All standards, controls and samples should be run in duplicate.
6. Samples containing IR [pY1158] protein extracted from cells should be diluted at least 1:10 with *Standard Diluent Buffer*. This dilution is necessary to reduce the matrix effect of the cell lysis buffer.
7. When pipetting reagents, maintain a consistent order of addition from well-to-well. This ensures equal incubation times for all wells.
8. Cover or cap all reagents when not in use.
9. **Do not mix or interchange different reagent lots from various kit lots.**
10. Do not use reagents after the kit expiration date.
11. Read absorbances within 2 hours of assay completion.
12. In-house controls should be run with every assay. If control values fall outside pre-established ranges, the accuracy of the assay is suspect.
13. All residual wash liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. *Never* insert absorbent paper directly into the wells.
14. Because *Stabilized Chromogen* is light sensitive, avoid prolonged exposure to light. Also avoid contact between *Stabilized Chromogen* and metal, or color may develop.

SAFETY

All blood components and biological materials should be handled as potentially hazardous. Follow universal precautions as established by the Centers for Disease Control and Prevention and by the Occupational Safety and Health Administration when handling and disposing of infectious agents.

DIRECTIONS FOR WASHING

Incomplete washing will adversely affect the test outcome. All washing must be performed with *Wash Buffer Concentrate (25X)* provided.

Washing can be performed manually as follows: completely aspirate the liquid from all wells by gently lowering an aspiration tip (aspiration device) into the bottom of each well. Take care not to scratch the inside of the well.

After aspiration, fill the wells with at least 0.4 mL of diluted wash solution. Let soak for 15 to 30 seconds, then aspirate the liquid. Repeat as directed under **ASSAY METHOD**. After the washing procedure, the plate is inverted and tapped dry on absorbent tissue.

Alternatively, the wash solution may be put into a squirt bottle. If a squirt bottle is used, flood the plate with wash buffer, completely filling all wells. After the washing procedure, the plate is inverted and tapped dry on absorbent tissue.

If using an automated washer, the operating instructions for washing equipment should be carefully followed. If your automated washer allows, 30 second soak cycles should be programmed into the wash cycle.

PROCEDURE FOR EXTRACTION OF PROTEINS FROM CELLS

A. Recommended Formulation of Cell Extraction Buffer:

10 mM Tris, pH 7.4

100 mM NaCl

1 mM EDTA

1 mM EGTA

1 mM NaF

20 mM $\text{Na}_4\text{P}_2\text{O}_7$

2 mM Na_3VO_4

1% Triton X-100

10% glycerol

0.1% SDS

0.5% deoxycholate

1 mM PMSF (stock is 0.3 M in DMSO)

Protease inhibitor cocktail (e.g., Sigma Cat. # P-2714) (reconstituted according to manufacturer's guideline). Add 250 μL per 5 mL Cell Extraction Buffer.

This buffer is stable for 2 to 3 weeks at 4°C or for up to 6 months when aliquoted (without protease inhibitors and PMSF added) and stored at -20°C. When stored frozen, the Cell Extraction Buffer should be thawed on ice. This buffer (minus protease inhibitors cocktail) can be obtained from Invitrogen under catalog # FNN0011. Important: add the protease inhibitors just before using. The stability of protease inhibitor supplemented Cell Extraction Buffer is 24 hours at 4°C. PMSF is very unstable and must be added prior to use, even if added previously.

B. Protocol for Cell Extraction

This protocol has been successfully applied to several cell lines. Researchers should optimize the cell extraction procedures for their own applications.

1. Collect cells in PBS by centrifugation (non-adherent) or scraping from culture flasks (adherent).
2. Wash cells twice with cold PBS.
3. Remove and discard the supernatant and collect the cell pellet. (At this point the cell pellet can be frozen at -80°C and lysed at a later date).
4. Lyse the cell pellet in Cell Extraction Buffer for 30 minutes, on ice, with vortexing at 10 minute intervals. The volume of Cell Extraction Buffer depends on the cell number in cell pellet and expression of IR [pY1158].
5. Transfer extract to microcentrifuge tubes and centrifuge at 13,000 rpm for 10 minutes at 4°C .
6. Aliquot the clear lysate to clean microfuge tubes. These samples are ready for assay. Lysates can be stored at -80°C . Avoid multiple freeze-thaws.

REAGENT PREPARATION AND STORAGE

A. Reconstitution and Dilution of IR [pY1158] Standard

Note: The *IR [pY1158] Standard* was prepared from autophosphorylated recombinant protein. One Unit of standard is equivalent to the amount of IR [pY1158] autophosphorylated from 0.6 ng of full-length (β -subunit) protein.

1. Reconstitute *IR [pY1158] Standard* with *Standard Diluent Buffer*. Refer to standard vial label for instructions. Swirl or mix gently and allow to sit for 10 minutes to ensure complete reconstitution. Label as 100 Units/mL IR [pY1158]. Use standard within 1 hour of reconstitution.
2. Add 0.25 mL of *Standard Diluent Buffer* to each of 6 tubes labeled 50, 25, 12.5, 6.25, 3.12 and 1.6 Units/mL IR [pY1158].
3. Make serial dilutions of the standard as described in the following dilution table. Mix thoroughly between steps.

B. Dilution of IR [pY1158] Standard

Standard:	Add:	Into:
100 Units/mL	Prepare as described in step 1	
50 Units/mL	0.25 mL of the 100 Units/mL std.	0.25 mL of the Diluent Buffer
25 Units/mL	0.25 mL of the 50 Units/mL std.	0.25 mL of the Diluent Buffer
12.5 Units/mL	0.25 mL of the 25 Units/mL std.	0.25 mL of the Diluent Buffer
6.25 Units/mL	0.25 mL of the 12.5 Units/mL std.	0.25 mL of the Diluent Buffer
3.12 Units/mL	0.25 mL of the 6.25 Units/mL std.	0.25 mL of the Diluent Buffer
1.6 Units/mL	0.25 mL of the 3.12 Units/mL std.	0.25 mL of the Diluent Buffer
0 Units/mL	0.25 mL of the Diluent Buffer	An empty tube

Discard all remaining reconstituted and diluted standards after completing assay. Return the *Standard Diluent Buffer* to the refrigerator.

C. Storage and Final Dilution of Anti-Rabbit IgG HRP (100X)

Please Note: The *Anti-Rabbit IgG HRP (100X)* is in 50% glycerol. This solution is viscous. To ensure accurate dilution, allow *Anti-Rabbit IgG HRP (100X)* to reach room temperature. Gently mix. Pipette *Anti-Rabbit IgG HRP (100X)* slowly. Remove excess concentrate solution from pipette tip by gently wiping with clean absorbent paper.

1. Dilute 10 μL of this 100X concentrated solution with 1 mL of *HRP Diluent* for each 8-well strip used in the assay. Label as Anti-Rabbit IgG HRP Working Solution.

For Example:

<u># of 8-Well Strips</u>	<u>Volume of Anti-Rabbit IgG HRP (100X)</u>	<u>Volume of Diluent</u>
2	20 μL solution	2 mL
4	40 μL solution	4 mL
6	60 μL solution	6 mL
8	80 μL solution	8 mL
10	100 μL solution	10 mL
12	120 μL solution	12 mL

2. Return the unused *Anti-Rabbit IgG HRP (100X)* to the refrigerator.

D. Dilution of Wash Buffer

Allow the *Wash Buffer Concentrate (25X)* to reach room temperature and mix to ensure that any precipitated salts have redissolved. Dilute 1 volume of the *Wash Buffer Concentrate (25X)* with 24 volumes of deionized water (e.g., 50 mL may be diluted up to 1.25 liters, 100 mL may be diluted up to 2.5 liters). Label as Working Wash Buffer.

Store both the concentrate and the Working Wash Buffer in the refrigerator. The diluted buffer should be used within 14 days.

ASSAY METHOD: PROCEDURE AND CALCULATIONS

Be sure to read the *Procedural Notes/Lab Quality Control* section before carrying out the assay.

Allow all reagents to reach room temperature before use. Gently mix all liquid reagents prior to use.

Note: A standard curve must be run with each assay.

1. Determine the number of 8-well strips needed for the assay. Insert these in the frame(s) for current use. (Re-bag extra strips and frame. Store these in the refrigerator for future use.)
2. Add 100 μL of the *Standard Diluent Buffer* to zero wells. Well(s) reserved for chromogen blank should be left empty.
3. Add 100 μL of standards to the appropriate microtiter wells. Samples prepared in Cell Extraction Buffer must be diluted 1:10 or greater in *Standard Diluent Buffer* (for example, 10 μL sample into 90 μL buffer). While a 1:10 sample dilution has been found to be satisfactory, higher dilutions such as 1:25 or 1:50 may be optimal for sample resolution. The dilution chosen should be optimized for

each investigation's system. Tap gently on side of plate to thoroughly mix. (See **REAGENT PREPARATION AND STORAGE**, Section B.)

4. Cover wells with *plate cover* and incubate for **2 hours at room temperature**.
5. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
6. Pipette 100 μL of *IR [pY1158] Detection Antibody* solution into each well except the chromogen blank(s). Tap gently on the side of the plate to mix.
7. Cover wells with *plate cover* and incubate for **1 hour at room temperature**.
8. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
9. Add 100 μL *Anti-Rabbit IgG HRP Working Solution* to each well except the chromogen blank(s). (Prepare the working dilution as described in **REAGENT PREPARATION AND STORAGE**, Section C.)
10. Cover wells with the *plate cover* and incubate for **30 minutes at room temperature**.
11. Thoroughly aspirate or decant solution from wells and discard the liquid. Wash wells 4 times. See **DIRECTIONS FOR WASHING**.
12. Add 100 μL of *Stabilized Chromogen* to each well. The liquid in the wells will begin to turn blue.
13. Incubate for **30 minutes at room temperature and in the dark**.
Please Note: Do not cover the plate with aluminum foil or

metalized mylar. The incubation time for chromogen substrate is often determined by the microtiter plate reader used. Many plate readers have the capacity to record a maximum optical density (O.D.) of 2.0. The O.D. values should be monitored and the substrate reaction stopped before the O.D. of the positive wells exceed the limits of the instrument. The O.D. values at 450 nm can only be read after the *Stop Solution* has been added to each well. If using a reader that records only to 2.0 O.D., stopping the assay after 20 to 25 minutes is suggested.

14. Add 100 μL of *Stop Solution* to each well. Tap side of plate gently to mix. The solution in the wells should change from blue to yellow.
15. Read the absorbance of each well at 450 nm having blanked the plate reader against a chromogen blank composed of 100 μL each of *Stabilized Chromogen* and *Stop Solution*. Read the plate within 2 hours after adding the *Stop Solution*.
16. Plot on graph paper the absorbance of the standards against the standard concentration. (Optimally, the background absorbance may be subtracted from *all* data points, including standards, unknowns and controls, prior to plotting.) Draw the best smooth curve through these points to construct the standard curve. If using curve fitting software, the four parameter algorithm provides the best curve fit.
17. Read the IR [pY1158] concentrations for unknown samples and controls from the standard curve plotted in step 16. **Multiply value(s) obtained for sample(s) by dilution factor to correct for the dilution in step 3.** (Samples still producing signals higher than the highest standard (100 Units/mL) should be further diluted in *Standard Diluent Buffer* and reanalyzed, multiplying the concentration found by the appropriate dilution factor.)

18. Values of IR [pY1158] should be normalized for IR (β -subunit) content by parallel measurement with the Invitrogen IR (β -subunit) ELISA Kit (Cat. # KHR9111).

TYPICAL DATA

The following data were obtained for the various standards over the range of 0 to 100 Units/mL IR [pY1158].

<u>IR [pY1158] Standard (Units/mL)</u>	<u>Optical Density (450 nm)</u>
0	0.179
1.6	0.220
3.12	0.256
6.25	0.384
12.5	0.577
25	0.938
50	1.516
100	2.524

LIMITATIONS OF THE PROCEDURE

Do not extrapolate the standard curve beyond the 100 Units/mL standard point; the dose-response is non-linear in this region and accuracy is difficult to obtain. Dilute samples >100 Units/mL with *Standard Diluent Buffer*; reanalyze these and multiply results by the appropriate dilution factor.

The influence of various extraction buffers has not been thoroughly investigated. The rate of degradation of native IR [pY1158] in various matrices has not been investigated. Although IR [pY1158] degradation in the Cell Extraction Buffer described in this protocol has not been seen to date, the possibility of this occurrence cannot be excluded.

For Research Use Only. CAUTION: Not for human or animal therapeutic or diagnostic use.


PERFORMANCE CHARACTERISTICS

SENSITIVITY

The analytical sensitivity of this assay is <0.8 Units/mL of IR [pY1158]. This was determined by adding two standard deviations to the mean O.D. obtained when the zero standard was assayed 30 times.

The sensitivity of this ELISA was compared to Western blotting using cell lysate with known quantities of IR [pY1158]. The data presented in Figure 1 show that the sensitivity of the ELISA is about 2x greater than that of Western blotting. The bands shown in the Western blot data were developed using rabbit anti-IR [pY1158] and alkaline phosphatase conjugated anti-rabbit IgG followed by chemiluminescent substrate and autoradiography.

Figure 1

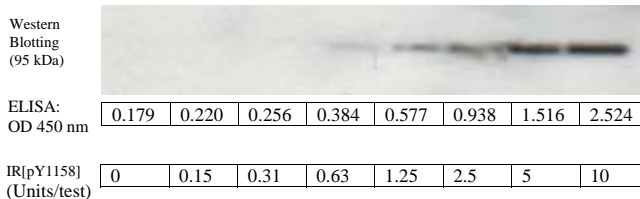

PRECISION

1. Intra-Assay Precision

Samples of known IR [pY1158] concentrations were assayed in replicates of 16 to determine precision within an assay.

	Sample 1	Sample 2	Sample 3
Mean (Units/mL)	66.36	18.58	7.65
SD	4.12	1.13	0.40
%CV	6.22	6.08	5.25

SD = Standard Deviation
CV = Coefficient of Variation

2. Inter-Assay Precision

Samples were assayed 48 times in multiple assays to determine precision between assays.

	Sample 1	Sample 2	Sample 3
Mean (Units/mL)	68.71	18.83	7.66
SD	5.33	1.60	0.48
%CV	7.76	8.48	6.29


SD = Standard Deviation
CV = Coefficient of Variation

RECOVERY

The recovery of IR [pY1158] added to 100 $\mu\text{g}/\text{mL}$ of a Jurkat cell lysate in Cell Extract Buffer (followed by 1:10 dilution in *Standard Diluent Buffer*) averaged 111%.

PARALLELISM

Natural IR [pY1158] from insulin-stimulated CHO-T was serially diluted in *Standard Diluent Buffer*. The optical density of each dilution was plotted against the IR [pY1158] standard curve. Parallelism was demonstrated by the figure below and indicated that the Standard accurately reflects IR [pY1158] content in samples.

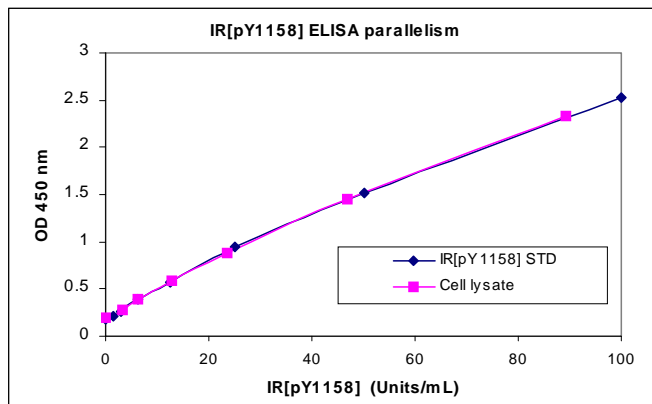

LINEARITY OF DILUTION


Extract Buffer was spiked with IR [pY1158] and serially diluted in *Standard Diluent Buffer* over the range of the assay. Linear regression analysis of samples versus the expected concentration yielded a correlation coefficient of 0.99.

Dilution	Cell Lysate		
	Measured (Units/mL)	Expected (Units/mL)	% Expected
Neat	89.1	-	-
1/2	46.8	44.6	105%
1/4	23.7	22.3	106%
1/8	12.8	11.1	114%
1/16	6.3	5.55	113%

SPECIFICITY

This IR [pY1158] ELISA kit is specific for measurement of IR that is phosphorylated at tyrosine 1158. IR of natural or recombinant (β -subunit) origin is reactive in this assay. This kit detects phosphorylated IR in insulin-stimulated CHO-T cells and does not detect non-phosphorylated IR in unstimulated cells, as shown in Figure 2.

Figure 2

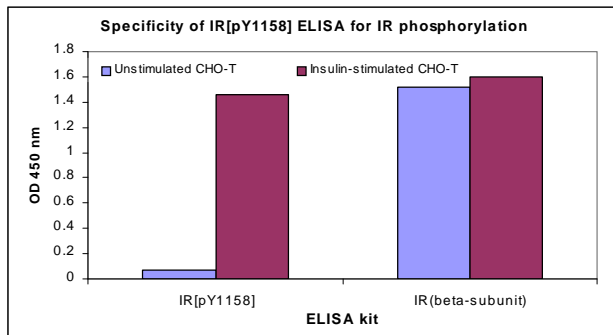

The specificity of this assay for IR phosphorylated at tyrosine 1158 was confirmed by peptide competition. Phosphorylated IR was quantitated in the assay as usual except that the detection antibody was preincubated with IR-derived peptides at a concentration of 0.01-1 μ g/mL. The data presented in Figure 3 show that only the

peptide corresponding to the region surrounding tyrosine 1158, containing the phospho-tyrosine, could block the ELISA signal. The peptides containing phosphorylated tyrosines at position 972, 1162/1163 did not block the signal.

Figure 3


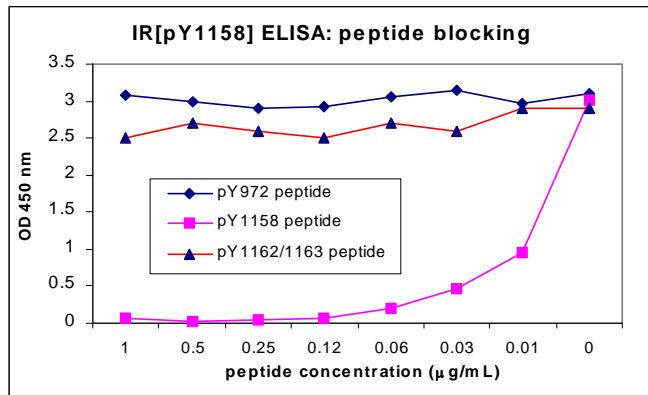


Figure 4 shows that IR [pY1158] phosphorylation in CHO-T cells is dependent on levels of insulin stimulation. Cells (approximately 90% confluent) were treated with insulin at varying concentrations (0-100 nM) for 10 minutes, lysed and quantitated in parallel for IR (β -subunit) and IR [pY1158] content. The amount of IR (β -subunit) remains constant, while the level of IR phosphorylated at tyrosine 1158 decreases with diminishing insulin dosage. Another IR phosphorylation site ELISA specific for [pYpY1162/1163] shows a similar pattern.

Figure 4




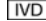


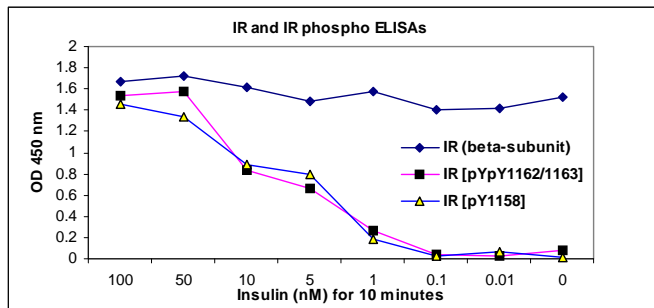








REFERENCES

1. Bevan, P., (2001) Insulin signaling. *J. Cell. Sci.* 114(Pt 8):1429-1430.
2. Skorey, K.L., et al. (2001) Development of a robust scintillation proximity assay for protein tyrosine phosphatase 1b using the catalytically inactive (c215s) mutant. *Anal. Biochem.* 291(2):269-278.
3. Ottensmeyer, F.P., et al. (2000) Mechanism of transmembrane signaling: insulin binding and the insulin receptor. *Biochemistry* 39(40):12103-12112.
4. Playford, M.P., et al. (2000) Insulin-like growth factor 1 regulates the location, stability, and transcriptional activity of beta-catenin. *Proc. Nat'l. Acad. Sci. USA* 97(22):12103–12108.
5. Virkamaki, A., et al. (1999) Protein-protein interaction in insulin signaling and the molecular mechanisms of insulin resistance. *J. Clin. Invest.* 103(7):931-943.
6. Wei, L., et al. (1995) Expression, characterization, and crystallization of the catalytic core of the human insulin receptor protein-tyrosine kinase domain. *J. Biol. Chem.* 270(14):8122-8130.
7. Ebina, Y., et al. (1985) The human insulin receptor cDNA: the structural basis for hormone-activated transmembrane signalling. *Cell* 40(4):747-758.
8. Ullrich, A., et al. (1985) Human insulin receptor and its relationship to the tyrosine kinase family of oncogenes. *Nature* 313(6005):756-761.

Important Licensing Information - These products may be covered by one or more Limited Use Label Licenses (see the Invitrogen Catalog or our website, www.invitrogen.com). By use of these products you accept the terms and conditions of all applicable Limited Use Label Licenses. Unless otherwise indicated, these products are for research use only and are not intended for human or animal diagnostic, therapeutic or commercial use.

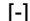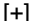Explanation of symbols

Symbol	Description	Symbol	Description
	Catalogue Number		Batch code
	Research Use Only		<i>In vitro</i> diagnostic medical device
	Use by		Temperature limitation
	Manufacturer		European Community authorised representative
	Without, does not contain		With, contains
	Protect from light		Consult accompanying documents
	Directs the user to consult instructions for use (IFU), accompanying the product.		

Copyright © Invitrogen Corporation. 22 February 2010

NOTES

Insulin Receptor [pY1158] Assay Summary

Incubate 100 μ L Standard or Cell Extract (>1:10)
for 2 hours at RT



aspirate and wash 4x

Incubate 100 μ L of Detection Antibody
for 1 hour at RT



aspirate and wash 4x

Incubate 100 μ L of HRP Anti-Rabbit Antibody
for 30 minutes at RT



aspirate and wash 4x

Incubate 100 μ L of Stabilized Chromogen
for 30 minutes at RT



Add 100 μ L of Stop Solution and read at 450 nm
Total time: 4 hours

